

USER GUIDE

Chromium Single Cell 3' Reagent Kits v3

with Feature Barcoding technology for
CRISPR Screening



FOR USE WITH

Chromium Single Cell 3' GEM, Library & Gel Bead Kit v3, 16 rxns PN-1000075

Chromium Single Cell 3' GEM, Library & Gel Bead Kit v3, 4 rxns PN-1000092

Chromium Single Cell 3' Feature Barcode Library Kit, 16 rxns PN-1000079

Chromium Single Cell B Chip Kit, 48 rxns PN-1000073

Chromium Single Cell B Chip Kit, 16 rxns PN-1000074

Chromium i7 Multiplex Kit, 96 rxns PN-120262

Notices

Document Number

CG000184 | Rev A

Legal Notices

© 2018 10X Genomics, Inc (10x Genomics). All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of 10x Genomics, is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any products described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. 10x Genomics provides no warranty and hereby disclaims any and all warranties as to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. A non-exhaustive list of 10x Genomics' marks, many of which are registered in the United States and other countries can be viewed at: www.10xgenomics.com/trademarks. 10x Genomics may refer to the products or services offered by other companies by their brand name or company name solely for clarity, and does not claim any rights in those third party marks or names. 10x Genomics products may be covered by one or more of the patents as indicated at: www.10xgenomics.com/patents. The use of products described herein is subject to 10x Genomics Terms and Conditions of Sale, available at www.10xgenomics.com/legal-notices, or such other terms that have been agreed to in writing between 10x Genomics and user. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

Instrument & Licensed Software Updates Warranties

Updates to existing Instruments and Licensed Software may be required to enable customers to use new or existing products. In the event of an Instrument failure resulting from an update, such failed Instrument will be replaced or repaired in accordance with the 10x Limited Warranty, Assurance Plan or service agreement, only if such Instrument is covered by any of the foregoing at the time of such failure. Instruments not covered under a current 10x Limited Warranty, Assurance Plan or service agreement will not be replaced or repaired.

Support

Email: support@10xgenomics.com

10x Genomics

7068 Koll Center Parkway

Suite 401

Pleasanton, CA 94566 USA

Table of Contents

Introduction	5
Chromium Single Cell 3' Reagent Kits v3	6
Chromium Accessories	10
Recommended Thermal Cyclers	10
Additional Kits, Reagents & Equipment	11
Protocol Steps & Timing	13
Stepwise Objectives	14
CRISPR Screening Overview	18
Tips & Best Practices	19
Step 1	25
GEM Generation & Barcoding	26
1.1 Prepare Master Mix	27
1.3 Run the Chromium Controller	30
1.4 Transfer GEMs	30
1.5 GEM-RT Incubation	31
Step 2	32
Post GEM-RT Cleanup & cDNA Amplification	33
2.1 Post GEM-RT Cleanup – Dynabeads	34
2.2 cDNA Amplification	36
2.3 cDNA Cleanup – SPRIselect	38
2.3A Pellet Cleanup	38
2.3B Transferred Supernatant Cleanup	38
2.4 Post cDNA Amplification QC & Quantification	39
Step 3	40
3' Gene Expression Library Construction	41
3.1 Fragmentation, End Repair & A-tailing	42
3.2 Post Fragmentation, End Repair & A-tailing Double Sided Size Selection – SPRIselect	43
3.3 Adaptor Ligation	44
3.4 Post Ligation Cleanup – SPRIselect	45
3.5 Sample Index PCR	46
3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect	47
3.7 Post Library Construction QC	48

Step 4	49
CRISPR Screening Library Construction	50
4.1 Guide RNA cDNA Cleanup – SPRIselect	51
4.2 Feature PCR	52
4.3 Post Feature PCR Cleanup – SPRIselect	53
4.4 Sample Index PCR	54
4.5 Post Sample Index PCR Double Sided Size Selection – SPRIselect	55
4.6 Post Library Construction QC	56
Sequencing	57
6.1 GEMs	61
6.2 Chromium Controller Errors	63
Appendix	64
Post Library Construction Quantification	65
Agilent TapeStation Traces	66
Compatible sgRNA Specifications	67
Oligonucleotide Sequences	68

Introduction

[Chromium Single Cell 3' Reagent Kits v3](#)

[Chromium Accessories](#)

[Recommended Thermal Cyclers](#)

[Additional Kits, Reagents & Equipment](#)

[Protocol Steps & Timing](#)

[Stepwise Objectives](#)

[CRISPR Screening Overview](#)

Chromium Single Cell 3' Reagent Kits v3

Chromium Single Cell 3' GEM, Library & Gel Bead Kit v3, 16 rxns PN-1000075

Chromium Single Cell 3' GEM Kit v3 16 rxns PN-1000077 (store at -20°C)

Chromium Single Cell 3' GEM Module

	#	PN
<input checked="" type="radio"/> RT Reagent	1	2000086
<input checked="" type="radio"/> RT Enzyme C	1	2000085
<input checked="" type="radio"/> Template Switch Oligo	1	3000228
<input type="radio"/> Reducing Agent B	1	2000087
<input type="radio"/> Cleanup Buffer	2	2000088
<input checked="" type="radio"/> cDNA Primers	1	2000089
<input type="radio"/> Amp Mix	1	2000047

10xGenomics.com

10x
GENOMICS

Chromium Single Cell 3' Library Kit v3 16 rxns PN-1000078 (store at -20°C)

Chromium Single Cell 3' Library Module

	#	PN
<input checked="" type="radio"/> Fragmentation Enzyme	1	2000090
<input checked="" type="radio"/> Fragmentation Buffer	1	2000091
<input checked="" type="radio"/> Ligation Buffer	1	2000092
<input checked="" type="radio"/> DNA Ligase	1	220110
<input checked="" type="radio"/> Adaptor Oligos	1	2000094
<input checked="" type="radio"/> SI Primer	1	2000095
<input type="radio"/> Amp Mix	1	2000047

10xGenomics.com

10x
GENOMICS

Chromium Single Cell 3' Gel Bead Kit v3, 16 rxns PN-1000076 (store at -80°C)

Chromium Single Cell 3' v3 Gel Beads

	#	PN
Single Cell 3' v3 Gel Beads	2	2000059

10xGenomics.com

10x
GENOMICS








Dynabeads™ MyOne™ SILANE PN-2000048 (store at 4°C)

	#	PN
Dynabeads MyOne SILANE	1	2000048

Chromium Single Cell 3' GEM, Library & Gel Bead Kit v3, 4 rxns PN-1000092

Chromium Single Cell 3' GEM Kit v3 4 rxns PN-1000094 (store at -20°C)

Chromium Single Cell 3' GEM Module







	#	PN
 RT Reagent	1	2000086
 RT Enzyme C	1	2000102
 Template Switch Oligo	1	3000228
 Reducing Agent B	1	2000087
 Cleanup Buffer	1	2000088
 cDNA Primers	1	2000089
 Amp Mix	1	2000103

10xGenomics.com

10x
GENOMICS

Chromium Single Cell 3' Library Kit v3 4 rxns PN-1000095 (store at -20°C)

Chromium Single Cell 3' Library Module

	#	PN
 Fragmentation Enzyme	1	2000104
 Fragmentation Buffer	1	2000091
 Ligation Buffer	1	2000092
 DNA Ligase	1	220131
 Adaptor Oligos	1	2000094
 SI Primer	1	2000095

10xGenomics.com

10x
GENOMICS

Chromium Single Cell 3' Gel Bead Kit v3, 4 rxns PN-1000093 (store at -80°C)

Chromium Single Cell 3' v3 Gel Beads

	#	PN
Single Cell 3' v3 Gel Beads (4 rxns)	1	2000059

10xGenomics.com


10x
GENOMICS

Dynabeads™ MyOne™ SILANE PN-2000048 (store at 4°C)

	#	PN
Dynabeads MyOne SILANE	1	2000048

Chromium Single Cell 3' Feature Barcode Library Kit, 16 rxns PN-1000079 (store at -20°C)


Chromium Single Cell 3' Feature Barcode Library Kit		
	#	PN
● Feature cDNA Primers 1	1	2000096
● Feature cDNA Primers 2	1	2000097
● Feature SI Primers 1	1	2000098
● Feature SI Primers 2	1	2000099
○ Amp Mix	2	2000047

10xGenomics.com 

Chromium Chip B Single Cell Kit, 48 rxns PN-1000073 (store at ambient temperature)

Chromium Partitioning Oil			Chromium Recovery Agent		
	#	PN		#	PN
<input checked="" type="radio"/> Partitioning Oil	6	220088	<input type="radio"/> Recovery Agent	6	220016


Chromium Chip B & Gaskets		
	#	PN
Chip B Single Cell	6	2000060
Gasket, 6-pack	1	370017

10xGenomics.com 

Chromium Chip B Single Cell Kit, 16 rxns PN-1000074 (store at ambient temperature)

Chromium Partitioning Oil			Chromium Recovery Agent		
	#	PN		#	PN
<input checked="" type="radio"/> Partitioning Oil	2	220088	<input type="radio"/> Recovery Agent	2	220016

Chromium Chip B & Gaskets		
	#	PN
Chip B Single Cell	2	2000060
Gasket, 2-pack	1	3000072

10xGenomics.com 

Chromium i7 Multiplex Kit, 96 rxns PN-120262 (store at -20°C)

Chromium i7 Multiplex Kit		
	#	PN
Chromium i7 Sample Index Plate	1	220103

Chromium Accessories

Product	Part Number (Kit)	Part Number (Item)
10x Vortex Adapter	120251	330002
10x Vortex Clip	120253	230002
10x Chip Holder	120252	330019
10x Magnetic Separator	120250	230003

Recommended Thermal Cyclers

Thermal cyclers used must support uniform heating of 100 μ l emulsion volumes.

Supplier	Description	Part Number
BioRad	C1000 Touch Thermal Cycler with 96-Deep Well Reaction Module	1851197
Eppendorf	MasterCycler Pro	North America 950030010 International 6321 000.019
Thermo Fisher Scientific	Veriti 96-Well Thermal Cycler	4375786

Additional Kits, Reagents & Equipment

The items in the table below have been validated by 10x Genomics and are highly recommended for the Chromium Single Cell 3' v3 protocols. Substituting materials may adversely affect system performance.







Supplier	Description	Part Number (US)
Plastics		
Eppendorf	PCR Tubes 0.2 ml 8-tube strips DNA LoBind Tubes, 1.5 ml DNA LoBind Tubes, 2.0 ml	951010022 022431021 022431048
USA Scientific	TempAssure PCR 8-tube strip	1402-4700
Thermo Fisher Scientific	MicroAmp 8-Tube Strip, 0.2 ml MicroAmp 8 -Cap Strip, clear	N8010580 N8010535
Rainin	Tips LTS 200UL Filter RT-L200FLR Tips LTS 1ML Filter RT-L1000FLR Tips LTS 20UL Filter RT-L10FLR	30389240 30389213 30389226
Kits & Reagents		
Thermo Fisher Scientific	Nuclease-free Water	AM9937
	Low TE Buffer (10 mM Tris-HCl pH 8.0, 0.1 mM EDTA)	12090-015
Millipore Sigma	Ethanol, Pure (200 Proof, anhydrous)	E7023-500ML
Beckman Coulter	SPRIselect Reagent Kit	B23318
Bio-Rad	10% Tween 20	1662404
Ricca Chemical Company	Glycerin (glycerol), 50% (v/v) Aqueous Solution	3290-32
Qiagen	Qiagen Buffer EB	19086
Equipment		
VWR	Vortex Mixer Divided Polystyrene Reservoirs	10153-838 41428-958
Eppendorf	Eppendorf ThermoMixer C Eppendorf SmartBlock 1.5 ml, Thermoblock for 24 reaction vessel (alternatively, use a temperature-controlled Heat Block)	5382000023 5360000038
Rainin	Pipet-Lite Multi Pipette L8-50XLS+ Pipet-Lite Multi Pipette L8-200XLS+ Pipet-Lite Multi Pipette L8-10XLS+ Pipet-Lite Multi Pipette L8-20XLS+ Pipet-Lite LTS Pipette L-2XLS+ Pipet-Lite LTS Pipette L-10XLS+ Pipet-Lite LTS Pipette L-20XLS+ Pipet-Lite LTS Pipette L-100XLS+ Pipet-Lite LTS Pipette L-200XLS+ Pipet-Lite LTS Pipette L-1000XLS+	17013804 17013805 17013802 17013803 17014393 17014388 17014392 17014384 17014391 17014382

Additional Kits, Reagents & Equipment

The items in the table below have been validated by 10x Genomics and are highly recommended for the Chromium Single Cell 3' v3 protocols. Substituting materials may adversely affect system performance.

Supplier	Description	Part Number (US)
Quantification & Quality Control		
Agilent	2100 Bioanalyzer Laptop Bundle	G2943CA
	High Sensitivity DNA Kit	5067-4626
	4200 TapeStation	G2991AA
	High Sensitivity D1000 ScreenTape/Reagents	5067-5592/ 5067-5593
	High Sensitivity D5000 ScreenTape/Reagents	5067-5584/ 5067-5585
		Choose Bioanalyzer, TapeStation, or Qubit based on availability & preference.
Thermo Fisher Scientific	Qubit 4.0 Fluorometer	Q33226
	Qubit dsDNA HS Assay Kit	Q32854
Advanced Analytical	Fragment Analyzer Automated CE System - 12 cap	FSv2-CE2F
	Fragment Analyzer Automated CE System - 48/96 cap	FSv2-CE10F
	High Sensitivity NGS Fragment Analysis Kit	DNF-474
KAPA Biosystems	KAPA Library Quantification Kit for Illumina Platforms	KK4824

Protocol Steps & Timing

Day	Steps	Timing	Stop & Store
2 h	Cell Preparation		
	Dependent on Cell Type	~1-1.5 h	
4 h	Step 1 – GEM Generation & Barcoding		
	1.1 Prepare Reaction Mix	20 min	
	1.2 Load Chromium Single Cell B Chip	10 min	
	1.3 Run the Chromium Controller	8.5 min	
	1.4 Transfer GEMs	3 min	
	1.5 GEM-RT Incubation	55 min	 4°C ≤72 h or -20°C ≤1 week
6 h	Step 2 – Post GEM-RT Cleanup & cDNA Amplification		
	2.1 Post GEM RT-Cleanup – Dynabead	45 min	
	2.2 cDNA Amplification	40 min	 4°C ≤72 h or -20°C ≤1 week
	2.3 cDNA Cleanup – SPRIselect		
	2.3A Pellet Cleanup	20 min	 4°C ≤72 h or -20°C ≤4 weeks
	2.3B Transferred Supernatant Cleanup	30 min	 4°C ≤72 h or -20°C ≤4 weeks
2.4 cDNA QC & Quantification	50 min		
8 h Plus	Step 3 – 3' Gene Expression Library Construction		
	3.1 Fragmentation, End Repair & A-tailing	50 min	
	3.2 Post Fragmentation, End Repair & A-tailing Double Sided Size Selection – SPRIselect	30 min	
	3.3 Adaptor Ligation	25 min	
	3.4 Post Ligation Cleanup – SPRIselect	20 min	
	3.5 Sample Index PCR	40 min	 4°C ≤72 h
	3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect	30 min	 4°C ≤72 h or -20°C long term
	3.7 Post Library Construction QC	50 min	
8 h Plus	Step 4 – CRISPR Screening Library Construction		
	4.1 Guide RNA cDNA Cleanup	20 min	
	4.2 Feature PCR	50 min	
	4.3 Post Feature PCR Cleanup – SPRIselect	20 min	
	4.4 Sample Index PCR	30 min	
	4.5 Post Sample Index PCR Size Selection – SPRIselect	20 min	 4°C ≤72 h or -20°C long term
	4.6 Post Library Construction QC	50 min	

Stepwise Objectives




The Chromium Single Cell Gene Expression Solution with Feature Barcoding technology upgrades short read sequencers to deliver a scalable microfluidic platform for assessing CRISPR-mediated perturbations of gene expression by profiling 500–10,000 individual cells per sample. GemCode Technology samples a pool of ~3,500,000 10x Barcodes to separately index each cell's transcriptome along with the CRISPR-mediated perturbations. It does so by partitioning thousands of cells into nanoliter-scale Gel Beads-in-emulsion (GEMs), where all generated cDNA (from poly-adenylated mRNAs and single-guide RNAs/sgRNAs) share a common 10x Barcode. Libraries are generated and sequenced from the cDNA and 10x Barcodes are used to associate individual reads back to the individual partitions.

This document outlines the protocol for generating Single Cell 3' Gene Expression and CRISPR Screening libraries from the same cells.

Single Cell 3' v3 Gel Beads


In addition to the poly(dT) primer that enables the production of barcoded, full-length cDNA from poly-adenylated mRNA, the Single Cell 3' v3 Gel Beads also include two additional primer sequences (Capture Sequence 1 and Capture Sequence 2), that enable capture and priming of Feature Barcoding technology compatible targets or analytes of interest.

The poly(dT) primers along with one of the capture sequence primers are used in this protocol for generating Single Cell 3' Gene Expression and CRISPR Screening libraries.

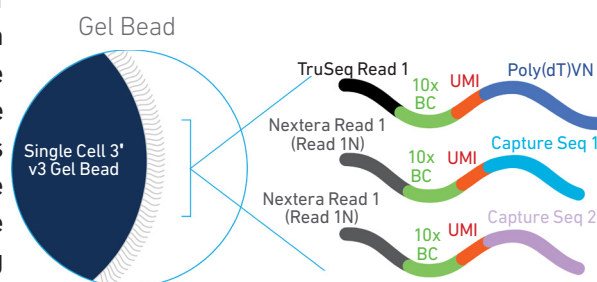

Step 1 GEM Generation & Barcoding

GEMs are generated by combining barcoded Single Cell 3' v3 Gel Beads, a Master Mix containing cells, and Partitioning Oil onto Chromium Chip B. To achieve single cell resolution, cells are delivered at a limiting dilution, such that the majority (~90-99%) of generated GEMs contain no cell, while the remainder largely contain a single cell.

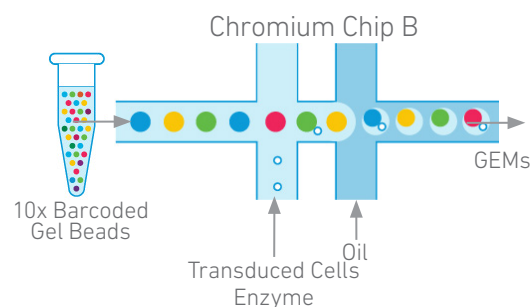

Step 1 GEM Generation & Barcoding

Immediately following GEM generation, the Gel Bead is dissolved releasing the three types of primers and any co-partitioned cell is lysed. The poly(dT) and the capture sequence primers in the gel bead are engaged simultaneously in two different reactions inside individual GEMs (primer with Capture Sequence 2 is not shown in the illustrated example).

A. Primers containing:

- an Illumina TruSeq Read 1 (read 1 sequencing primer)
- 16 nt 10x Barcode
- 12 nt unique molecular identifier (UMI)
- 30 nt poly(dT) sequence

are mixed with the cell lysate and a Master Mix containing reverse transcription (RT) reagents. Incubation of the GEMs produces barcoded, full-length cDNA from poly-adenylated mRNA.


B. In the same partition, primers containing:

- an Illumina Nextera Read 1 (Read 1N; read 1 sequencing primer)
- 16 nt 10x Barcode
- 12 nt unique molecular identifier (UMI)
- Capture Sequence 1

are mixed with the cell lysate and the Master Mix containing RT reagents. Incubation of the GEMs produces barcoded, cDNA from the sgRNA.

[See Appendix for Compatible sgRNA Specifications](#)

Inside individual GEMs

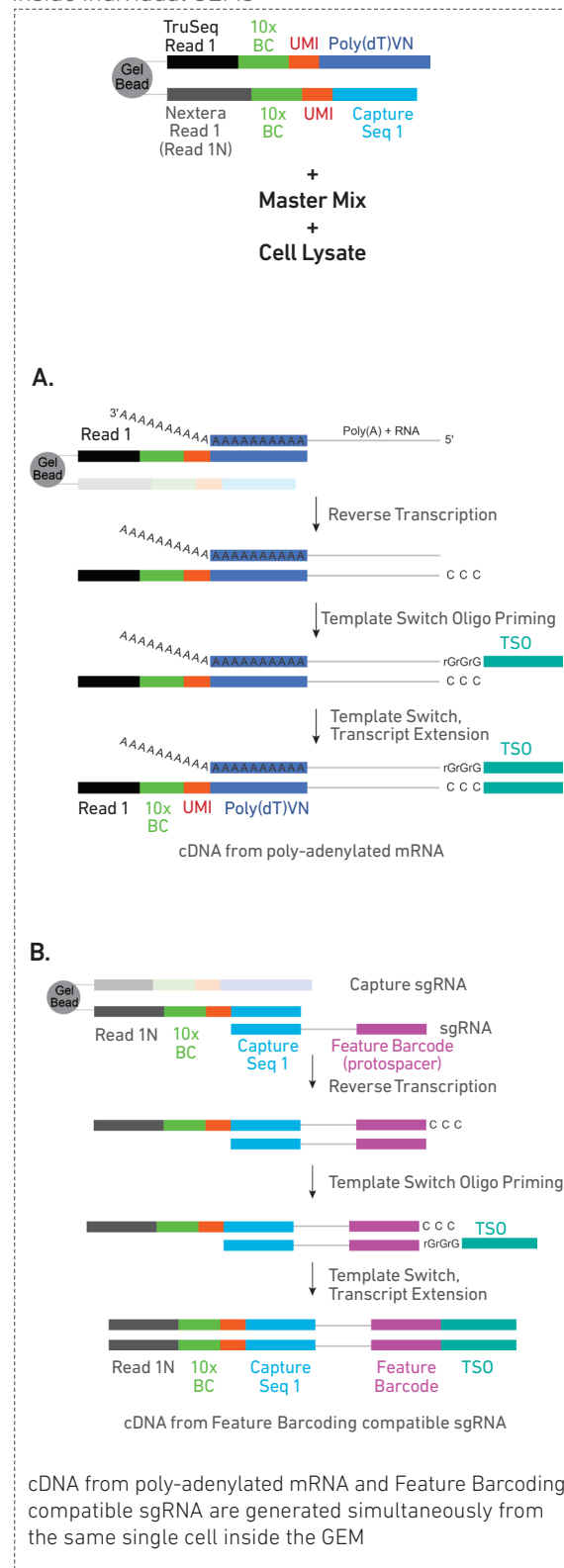


Step 2 Post GEM-RT Cleanup & cDNA Amplification



After incubation, GEMs are broken and pooled fractions are recovered. Silane magnetic beads are used to purify the cell barcoded products from the post GEM-RT reaction mixture, which includes leftover biochemical reagents and primers. The cell barcoded cDNA molecules are amplified via PCR to generate sufficient mass for library constructions. Size selection is used to separate the amplified cDNA molecules for 3' Gene Expression and CRISPR Screening library construction.

Pooled cDNA amplification

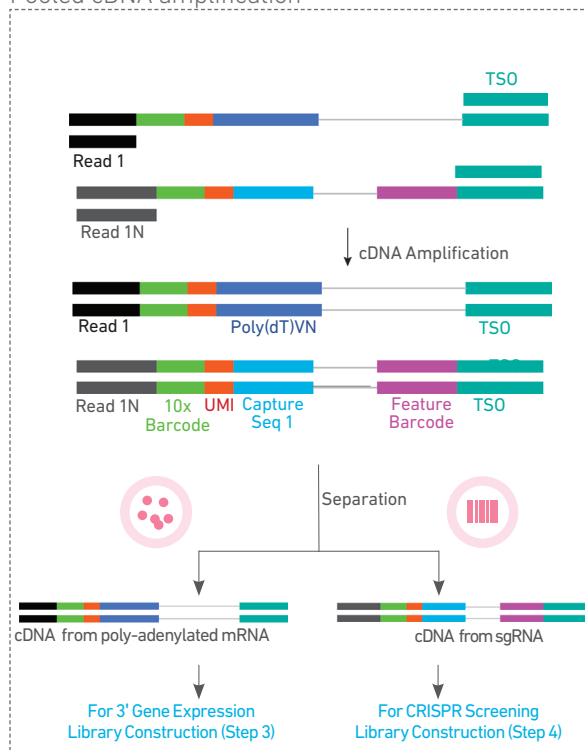


Step 3 3' Gene Expression Library Construction



Full-length cDNA from poly-adenylated mRNA is used to generate 3' Gene Expression libraries. Enzymatic fragmentation and size selection are used to optimize the cDNA amplicon size. TruSeq Read 1 (read 1 primer sequence) is added to the molecules during GEM incubation. P5, P7, a sample index, and TruSeq Read 2 (read 2 primer sequence) are added via End Repair, A-tailing, Adaptor Ligation, and PCR. The final libraries contain the P5 and P7 primers used in Illumina bridge amplification.

Pooled amplified cDNA processed in bulk

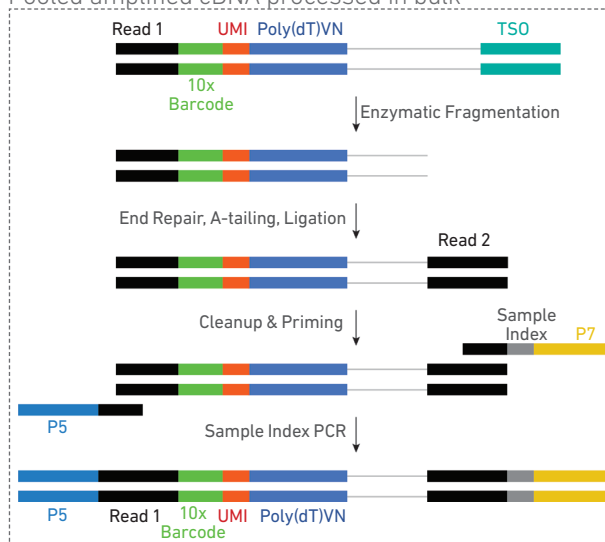


Step 4 CRISPR Screening Library Construction



Amplified cDNA from sgRNA molecules is used to generate CRISPR Screening libraries. P5, P7, a sample index, and TruSeq Read 2 (read 2 primer sequence) are added via PCR. The final libraries contain the P5 and P7 primers used in Illumina bridge amplification.

Pooled amplified DNA processed in bulk

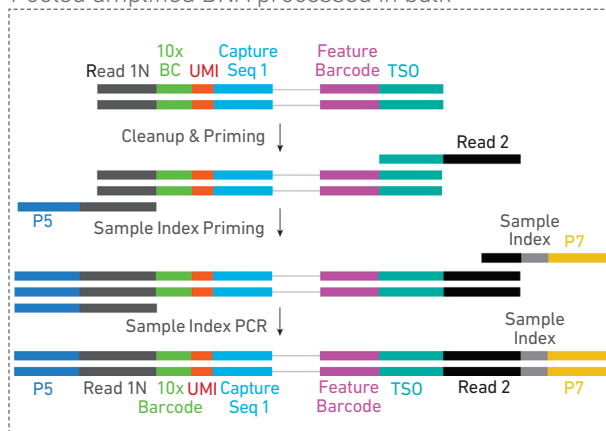


Step 5 Sequencing


The Single Cell 3' libraries comprise standard Illumina paired-end constructs which begin and end with P5 and P7. The 16 bp 10x Barcode and 12 bp UMI are encoded in Read 1, while Read 2 is used to sequence the cDNA fragment in 3' Gene Expression libraries and the Feature Barcode (sgRNA protospacer) in the CRISPR Screening libraries. Sample index sequences are incorporated as the i7 index read. Standard Illumina sequencing primer sites TruSeq Read 1 and TruSeq Read 2 in the 3' Gene Expression libraries and Nextera Read 1 and Truseq Read 2 in the CRISPR Screening libraries are used in paired-end sequencing.

Illumina sequencer compatibility, sample indices, library loading and pooling, recommended read depths & run parameters for sequencing are summarized in step 5.

Chromium Single Cell 3' Gene Expression Library

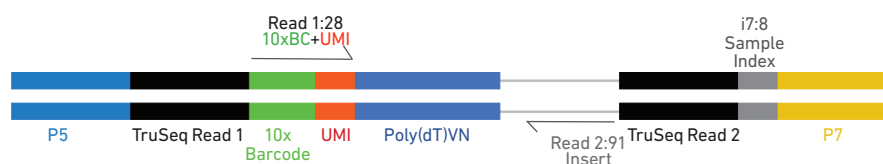

Chromium Single Cell 3' CRISPR Screening Library

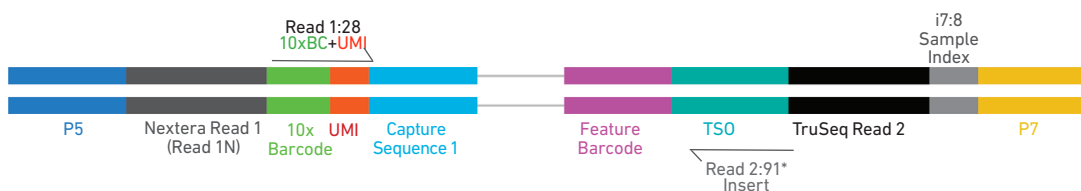


*Minimum required Read 2 length for CRISPR Screening libraries is 70 bp

[See Appendix for Oligonucleotide Sequences](#)

CRISPR Screening Overview


The Chromium Single Cell Gene Expression Solution with Feature Barcoding technology provides a high-throughput and scalable approach to obtain gene expression profiles along with perturbation phenotypes via direct capture of poly-adenylated mRNAs and single-guide RNAs (sgRNAs) from the same single cell (see [Stepwise Objectives](#)). For compatibility with Feature Barcoding technology, sgRNAs should be engineered containing either Capture Sequence 1 or Capture Sequence 2. Two possible locations for integrating the capture sequence in the sgRNA include (1) within the sgRNA hairpin structure, or (2) immediately before the sgRNA termination signal, elongating the 3'-end of the sgRNA. However, alternate sgRNA integration locations for either of the two capture sequences may be possible depending on the specific application, type of construct used etc.

Capture Sequence Integration in sgRNA

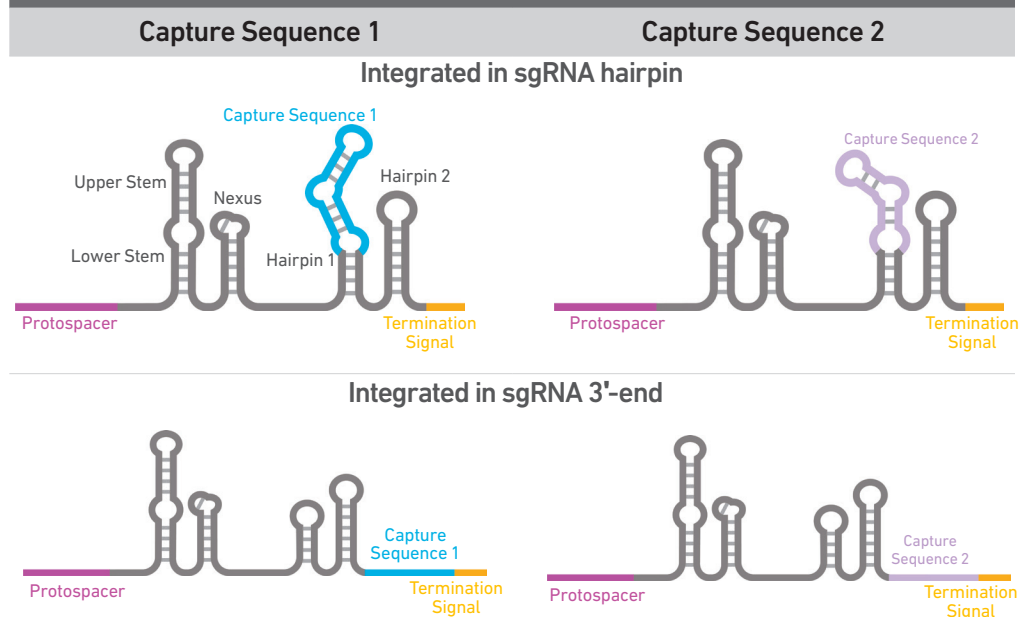

Direct Capture of sgRNA

Only Capture Sequence 1 is illustrated in the example

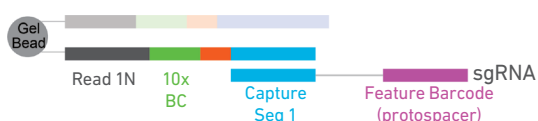

Performing sgRNA QC by qPCR, NGS, or other methods is recommended prior to proceeding with the Single Cell Gene Expression and CRISPR Screening Solution.

[See Appendix for Compatible sgRNA Specifications](#)

Tips & Best Practices



Icons



Tips & Best Practices section includes additional guidance



Signifies critical step requiring accurate execution



Troubleshooting section includes additional guidance

Version Specific Update



Indicates version specific updates in a particular protocol step to inform users who have used a previous version of the product. The updates may be in volume, temperature, calculation instructions etc.

Emulsion-safe Plastics

- Use validated emulsion-safe plastic consumables when handling GEMs as some plastics can destabilize GEMs.

Cell Concentration

- Recommended starting point is to load ~1,600 cells per reaction, resulting in recovery of ~1,000 cells, and a multiplet rate of ~0.8%. The optimal input cell concentration is 700-1,200 cells/ μ l.
- The presence of dead cells in the suspension may also reduce the recovery rate. Consult the 10x Genomics Single Cell Protocols Cell Preparation Guide and the Guidelines for Optimal Sample Preparation flowchart (Documents CG00053 and CG000126 respectively) for more information on preparing cells.

Multiplet Rate (%)	# of Cells Loaded	# of Cells Recovered
~0.4%	~800	~500
~0.8%	~1,600	~1,000
~1.6%	~3,200	~2,000
~2.3%	~4,800	~3,000
~3.1%	~6,400	~4,000
~3.9%	~8,000	~5,000
~4.6%	~9,600	~6,000
~5.4%	~11,200	~7,000
~6.1%	~12,800	~8,000
~6.9%	~14,400	~9,000
~7.6%	~16,000	~10,000

General Reagent Handling

- Fully thaw and thoroughly mix reagents before use.
- Keep all enzymes and Master Mixes on ice during setup and use. Promptly move reagents back to the recommended storage.
- Calculate reagent volumes with 10% excess of 1 reaction values.
- Cover Partitioning Oil tubes and reservoirs to minimize evaporation.
- If using multiple chips, use separate reagent reservoirs for each chip during loading.
- Thoroughly mix samples with the beads during bead-based cleanup steps.

50% Glycerol Solution

- Purchase 50% glycerol solution from Ricca Chemical Company, Glycerin (glycerol), 50% (v/v) Aqueous Solution, PN-3290-32. OR
- Prepare 50% glycerol solution:
 - i. Mix an equal volume of water and 99% Glycerol, Molecular Biology Grade.
 - ii. Filter through a 0.2 μm filter.
 - iii. Store at -20°C in 1-ml LoBind tubes. 50% glycerol solution should be equilibrated to room temperature before use.

Pipette Calibration


- Follow manufacturer's calibration and maintenance schedules.
- Pipette accuracy is particularly important when using SPRIselect reagents.

Chromium Chip Handling

- Minimize exposure of reagents, chips, and gaskets to sources of particles and fibers, laboratory wipes, frequently opened flip-cap tubes, clothing that sheds fibers, and dusty surfaces.
- Execute steps without pause or delay, unless indicated. When multiple chips are to be used, load, run, and collect the content from one chip before loading the next.
- Fill all unused input wells in rows labeled 1, 2, and 3 on a chip with an appropriate volume of 50% glycerol solution before loading the used wells. DO NOT add glycerol to the Recovery Wells.
- Avoid contacting the bottom surface of the chip with gloved hands and other surfaces. Frictional charging can lead to inadequate priming of the channels, potentially leading to either clogs or wetting failures.
- Minimize the distance that a loaded chip is moved to reach the Chromium Controller.
- Keep the chip horizontal to prevent wetting the gasket with oil, which depletes the input volume and may adversely affect the quality of the resulting emulsion.


10x Chip Holders

- 10x Chip Holders encase Chromium Chips.
- The holder lid flips over to become a stand, holding the chip at 45 degrees for optimal recovery well content removal.
- Squeeze the black sliders on the back side of the holder together to unlock the lid and return the holder to a flat position.

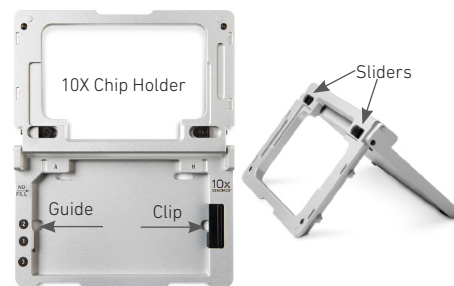

Chromium Chip & Holder Assembly

- Align notch on the chip (upper left corner) and the holder.
- Insert the left-hand side of the chip under the guide. Depress the right-hand side of the chip until the spring-loaded clip engages.
- Close the lid before dispensing reagents into the wells.

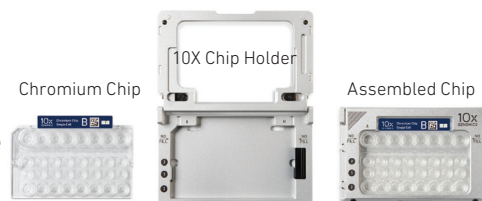

Chromium Chip Loading


- Place the assembled chip and holder flat on the bench with the lid closed.
- Dispense at the bottom of the wells without introducing bubbles.
- When dispensing Gel Beads into the chip, wait for the remainder to drain into the bottom of the pipette tips and dispense again to ensure complete transfer.
- Refer to [Load Chromium Chip B](#) for specific instructions.



Gel Bead Handling



- Use one tube of Gel Beads per sample. DO NOT puncture the foil seals of tubes not used at the time.
- Equilibrate the Gel Beads strip to room temperature before use.
- Store unused Gel Beads at -80°C and avoid more than 12 freeze-thaw cycles. DO NOT store Gel Beads at -20°C .
- Attach a 10x Vortex Adapter to the top of standard laboratory vortexers to vortex the Gel Bead strips.
- After vortexing, remove the Gel Bead strip from the adapter. Flick the Gel Bead strip in a sharp, downward motion maximize Gel Bead recovery. Confirm there are no bubbles at the bottom of the tubes.
- If the required volume of beads cannot be recovered, place the pipette tips against the sidewalls and slowly dispense the Gel Beads back into the tubes. DO NOT introduce bubbles into the tubes and verify that the pipette tips contain no leftover Gel Beads. Withdraw the full volume of beads again by pipetting slowly.

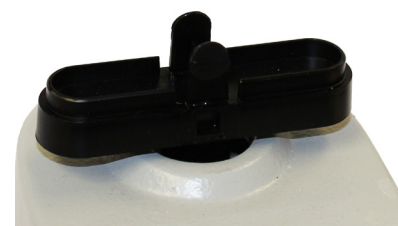

10x Gasket Attachment

- After reagents are loaded, attach the gasket by holding the tongue (curved end, to the right) and hook it on the left-hand tabs of the holder. Gently pull the gasket toward the right and hook it on the two right-hand tabs.
- DO NOT touch the smooth side of the gasket. DO NOT press down on the top of the gasket after attachment.
- Keep the assembly horizontal to avoid wetting the gasket with Partitioning Oil.



10x Magnetic Separator


- Offers two positions of the magnets (high and low) relative to a tube, depending on its orientation. Flip the magnetic separator over to switch between high (magnet•High) or low (magnet•Low) positions.
- If using MicroAmp 8-Tube Strips, use the high position (magnet•High) only throughout the protocol.




Magnetic Bead Cleanup Steps

- During magnetic bead based cleanup steps that specify waiting “until the solution clears”, visually confirm clearing of solution before proceeding to the next step. See adjacent panel for an example.
- The time needed for the solution to clear may vary based on specific step, reagents, volume of reagents etc.


Magnetic beads mixed with reagent

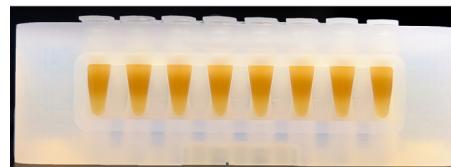


Separation in progress

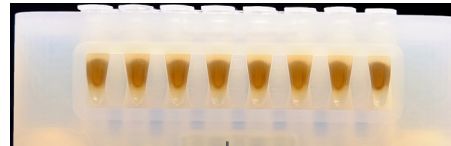

Separation continuing

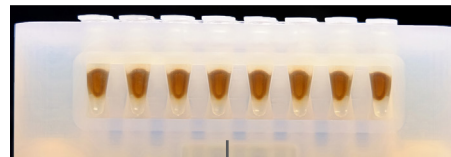

Separation complete; solution is clear

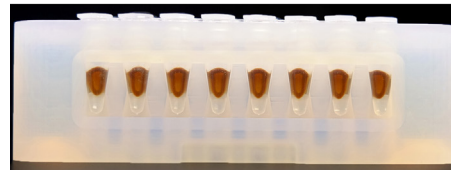

SPRIselect Cleanup & Size Selection


- After aspirating the desired volume of SPRIselect reagent, examine the pipette tips before dispensing to ensure the correct volume is transferred.
- Pipette mix thoroughly as insufficient mixing of sample and SPRIselect reagent will lead to inconsistent results.
- Use fresh preparations of 80% Ethanol.

Tutorial — SPRIselect Reagent:DNA Sample Ratios

SPRI beads selectively bind DNA according to the ratio of SPRIselect reagent (beads).

$$\text{Example: Ratio} = \frac{\text{Volume of SPRIselect reagent added to the sample}}{\text{Volume of DNA sample}} = \frac{50 \mu\text{l}}{100 \mu\text{l}} = 0.5\text{X}$$

Schematic of Double Sided Size Selection

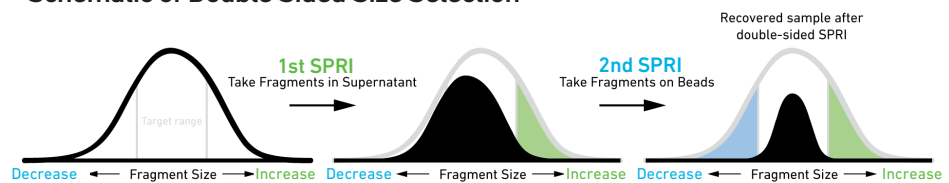

After the first SPRI, supernatant is transferred for a second SPRI while larger fragments are discarded (green). After the second SPRI, fragments on beads are eluted and kept while smaller fragments are discarded (blue). Final sample has a tight fragment size distribution with reduced overall amount (black).

Tutorial — Double Sided Size Selection

Step a – First SPRIselect: Add 50 μl SPRIselect reagent to 100 μl sample (0.5X).

$$\text{Ratio} = \frac{\text{Volume of SPRIselect reagent added to the sample}}{\text{Volume of DNA sample}} = \frac{50 \mu\text{l}}{100 \mu\text{l}} = 0.5\text{X}$$

Step b – Second SPRIselect: Add 30 μl SPRIselect reagent to supernatant from step a (0.8X).

$$\text{Ratio} = \frac{\text{Total Volume of SPRIselect reagent added to the sample (step a + b)}}{\text{Original Volume of DNA sample}} = \frac{50 \mu\text{l} + 30 \mu\text{l}}{100 \mu\text{l}} = 0.8\text{X}$$

Enzymatic Fragmentation

- Ensure enzymatic fragmentation reactions are prepared on ice and then loaded into a thermal cycler pre-cooled to 4°C prior to initiating the Fragmentation, End Repair, and A-tailing incubation steps.

Sample Indices in Sample Index PCR

- Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run.
- Each well in the i7 Sample Index plate contains a unique mix of 4 oligos.
- The sample indexes can therefore be used in any combination.
- Each sample index set is base-balanced to avoid monochromatic signal issues when it is the sole sample loaded on an Illumina sequencer.

Step 1


GEM Generation & Barcoding

- 1.1 Prepare Single Cell Master Mix
- 1.2 Load Chromium Chip B
- 1.3 Run the Chromium Controller
- 1.4 Transfer GEMs
- 1.5 GEM-RT Incubation

1

1.0
GEM Generation & Barcoding

VERSION SPECIFIC

GET STARTED!				
Action	Item	10x PN	Preparation & Handling	Storage
Equilibrate to Room Temperature	Chromium Single Cell 3' v3 Gel Beads	2000059	Equilibrate to room temperature 30 min before loading the chip.	-80°C
	<input checked="" type="radio"/> RT Reagent	2000086	Vortex, verify no precipitate, centrifuge briefly.	-20°C
	<input checked="" type="radio"/> Template Switch Oligo	3000228	Centrifuge briefly, resuspend in 80 µl Low TE Buffer. Vortex 15 sec at maximum speed, centrifuge briefly, leave at room temperature for ≥ 30 min. After resuspension, store at -80°C.	-20°C
	<input type="radio"/> Reducing Agent B	2000087	Vortex, verify no precipitate, centrifuge briefly.	-20°C
Place on Ice	<input checked="" type="radio"/> RT Enzyme C	2000085/ 2000102	Centrifuge briefly before adding to the mix.	-20°C
Feature Barcoding technology compatible Cell Suspension				
 Consult Technical Note Guide RNA Specifications Compatible with Feature Barcoding technology for CRISPR Screening (CG000197)				
Obtain	<input type="radio"/> Partitioning Oil	220088	-	Ambient
	Chromium Chip B Single Cell	2000060	-	Ambient
	10x Gasket	370017/ 3000072	See Tips & Best Practices.	Ambient
	10x Chip Holder	330019	See Tips & Best Practices.	Ambient
	10x Vortex Adapter	330002	See Tips & Best Practices.	Ambient
	50% glycerol solution If using <8 reactions	-	See Tips & Best Practices.	-



Firmware Version 3.16 or higher is required in the Chromium Controller or the Chromium Single Cell Controller used for Single Cell 3' protocols.

1.1 Prepare Master Mix

a. Prepare Master Mix on ice. Pipette mix 15x and centrifuge briefly.

Master Mix <i>Add reagents in the order listed</i>	PN	1X (μl)	4X + 10% (μl)	8X + 10% (μl)
● RT Reagent	2000086	20.0	88.0	176.0
● Template Switch Oligo	3000228	3.1	13.9	27.7
○ Reducing Agent B	2000087	2.0	8.7	17.3
● RT Enzyme C	2000085/ 2000102	8.3	36.6	73.1
Total	-	33.4	147.1	294.2

b. Add 33.4 μl Master Mix into each tube of a PCR 8-tube strip on ice.

Cell Suspension Volume Calculator Table (for step 1.2)

Volume of Transduced Cell Suspension Stock per reaction (µl) | Volume of Nuclease-free Water per reaction (µl)

Cell Stock Concentration (Cells/µl)	Targeted Cell Recovery											
	500	1000	2000	3000	4000	5000	6000	7000	8000	9000	10000	
100	8.0 38.6	16.0 30.6	32.0 14.6	n/a	n/a	n/a	n/a	n/a	n/a	n/a	n/a	n/a
200	4.0 42.6	8.0 38.6	16.0 30.6	24.0 22.6	32.0 14.6	40.0 6.6	n/a	n/a	n/a	n/a	n/a	n/a
300	2.7 43.9	5.3 41.3	10.7 35.9	16.0 30.6	21.3 25.3	26.7 19.9	32.0 14.6	37.3 9.3	42.7 3.9	n/a	n/a	n/a
400	2.0 44.6	4.0 42.6	8.0 38.6	12.0 34.6	16.0 30.6	20.0 26.6	24.0 22.6	28.0 18.6	32.0 14.6	36.0 10.6	40.0 6.6	40.0
500	1.6 45.0	3.2 43.4	6.4 40.2	9.6 37.0	12.8 33.8	16.0 30.6	19.2 27.4	22.4 24.2	25.6 21.0	28.8 17.8	32.0 14.6	32.0
600	1.3 45.3	2.7 43.9	5.3 41.3	8.0 38.6	10.7 35.9	13.3 33.3	16.0 30.6	18.7 27.9	21.3 25.3	24.0 22.6	26.7 19.9	26.7
700	1.1 45.5	2.3 44.3	4.6 42.0	6.9 39.7	9.1 37.5	11.4 35.2	13.7 32.9	16.0 30.6	18.3 28.3	20.6 26.0	22.9 23.7	22.9
800	1.0 45.6	2.0 44.6	4.0 42.6	6.0 40.6	8.0 38.6	10.0 36.6	12.0 34.6	14.0 32.6	16.0 30.6	18.0 28.6	20.0 26.6	20.0
900	0.9 45.7	1.8 44.8	3.6 43.0	5.3 41.3	7.1 39.5	8.9 37.7	10.7 35.9	12.4 34.2	14.2 32.4	16.0 30.6	17.8 28.8	17.8
1000	0.8 45.8	1.6 45.0	3.2 43.4	4.8 41.8	6.4 40.2	8.0 38.6	9.6 37.0	11.2 35.4	12.8 33.8	14.4 32.2	16.0 30.6	16.0
1100	0.7 45.9	1.5 45.1	2.9 43.7	4.4 42.2	5.8 40.8	7.3 39.3	8.7 37.9	10.2 36.4	11.6 35.0	13.1 33.5	14.5 32.1	14.5
1200	0.7 45.9	1.3 45.3	2.7 43.9	4.0 42.6	5.3 41.3	6.7 39.9	8.0 38.6	9.3 37.3	10.7 35.9	12.0 34.6	13.3 33.3	13.3
1300	0.6 46.0	1.2 45.4	2.5 44.1	3.7 42.9	4.9 41.7	6.2 40.4	7.4 39.2	8.6 38.0	9.8 36.8	11.1 35.5	12.3 34.3	12.3
1400	0.6 46.0	1.1 45.5	2.3 44.3	3.4 43.2	4.6 42.0	5.7 40.9	6.9 39.7	8.0 38.6	9.1 37.5	10.3 36.3	11.4 35.2	11.4
1500	0.5 46.1	1.1 45.5	2.1 44.5	3.2 43.4	4.3 42.3	5.3 41.3	6.4 40.2	7.5 39.1	8.5 38.1	9.6 37.0	10.7 35.9	10.7
1600	0.5 46.1	1.0 45.6	2.0 44.6	3.0 43.6	4.0 42.6	5.0 41.6	6.0 40.6	7.0 39.6	8.0 38.6	9.0 37.6	10.0 36.6	10.0
1700	0.5 46.1	0.9 45.7	1.9 44.7	2.8 43.8	3.8 42.8	4.7 41.9	5.6 41.0	6.6 40.0	7.5 39.1	8.5 38.1	9.4 37.2	9.4
1800	0.4 46.2	0.9 45.7	1.8 44.8	2.7 43.9	3.6 43.0	4.4 42.2	5.3 41.3	6.2 40.4	7.1 39.5	8.0 38.6	8.9 37.7	8.9
1900	0.4 46.2	0.8 45.8	1.7 44.9	2.5 44.1	3.4 43.2	4.2 42.4	5.1 41.5	5.9 40.7	6.7 39.9	7.6 39.0	8.4 38.2	8.4
2000	0.4 46.2	0.8 45.8	1.6 45.0	2.4 44.2	3.2 43.4	4.0 42.6	4.8 41.8	5.6 41.0	6.4 40.2	7.2 39.4	8.0 38.6	8.0

Grey boxes: Volumes that would exceed the allowable water volume in each reaction

Yellow boxes: Indicate a low transfer volume that may result in higher cell load variability

Blue boxes: Optimal range of cell stock concentration to maximize the likelihood of achieving the desired cell recovery target

1.2 Load Chromium Chip B

See Tips & Best Practices for chip handling instructions. When loading the chip, raising and depressing the pipette plunger should each take ~5 sec. When dispensing, raise the pipette tips at the same rate as the liquid is rising, keeping the tips slightly submerged.

TIPS

a. **Assemble Chromium Chip B in a 10x Chip Holder.** See Tips & Best Practices.

VERSION SPECIFIC

b. **Dispense 50% Glycerol Solution into Unused Chip Wells** (if < 8 samples per chip)
 i. 75 μ l to unused wells in row labeled 1. DO NOT add 50% glycerol solution to the top row of Recovery Wells. DO NOT use any substitute for 50% glycerol solution.
 ii. 40 μ l to unused wells in row labeled 2.
 iii. 280 μ l to unused wells in row labeled 3.


c. **Prepare Master Mix + Cell Suspension**

Refer to the Cell Suspension Volume Calculator Table and add the appropriate volume of nuclease-free water and corresponding volume of transduced single cell suspension to Master Mix for a total of 80 μ l in each tube. Gently pipette mix the single cell suspension before adding to the Master Mix.

VERSION SPECIFIC

d. **Load Row Labeled 1**

Gently pipette mix the Master Mix + Cell Suspension and using the same pipette tip, dispense 75 μ l Master Mix + Cell Suspension into the bottom center of each well in row labeled 1 without introducing bubbles.

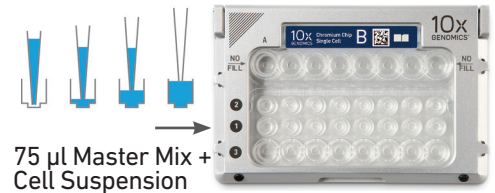


e. **Prepare Gel Beads**

Snap the Gel Bead strip into a 10x Vortex Adapter. Vortex 30 sec. Remove the Gel Bead strip and flick in a sharp, downward motion to ensure maximum recovery. Confirm there are no bubbles at the bottom of the tubes and liquid levels look even.

f. **Load Row Labeled 2**


Puncture the foil seal of the Gel Bead tubes. Slowly aspirate 40 μ l Gel Beads. Dispense into the wells in row labeled 2 without introducing bubbles.

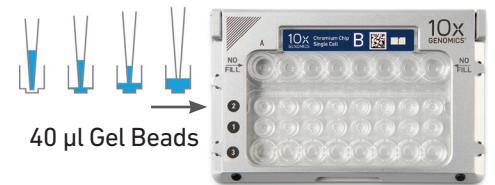

VERSION SPECIFIC

g. **Load Row Labeled 3**

Dispense 280 μ l Partitioning Oil into the wells in row labeled 3 by pipetting two aliquots of 140 μ l from a reagent reservoir. Failure to add Partitioning Oil can damage the Chromium Controller.

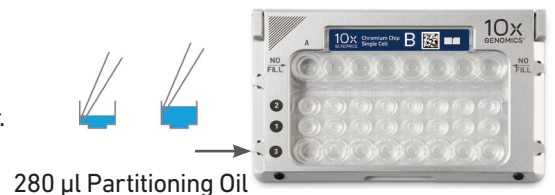

h. **Attach 10x Gasket**

Align the notch with the top left-hand corner. Ensure the gasket holes are aligned with the wells. Avoid touching the smooth gasket surface. DO NOT press down on the gasket.

Keep horizontal to avoid wetting the gasket.



1.3 Run the Chromium Controller

- Press the eject button on the Controller to eject the tray.
- Place the assembled chip with the gasket in the tray. Press the button to retract the tray.
- Confirm the Chromium Single Cell B program on screen. Press the play button.
- At completion of the run (~8.5 min), the Controller will chime. **Immediately** proceed to the next step.



1.4 Transfer GEMs

- Place a tube strip on ice.
- Press the eject button of the Controller and remove the chip.
- Discard the gasket. Open the chip holder. Fold the lid back until it clicks to expose the wells at 45 degrees. Ensure that the partitioning oil from the wells does not spill when exposing the wells.
- Check the volume in rows 1-3. Abnormally high volume in any well indicates a clog.
- Slowly aspirate **100 μ l** GEMs from the lowest points of the Recovery Wells in the top row without creating a seal between the tips and the bottom of the wells.
- Withdraw pipette tips from the wells. GEMs should appear opaque and uniform across all channels. Excess Partitioning Oil (clear) in the pipette tips indicates a potential clog.
- Over the course of **~20 sec**, dispense GEMs into the tube strip on ice with the pipette tips against the sidewalls of the tubes.
- If multiple chips are run back-to-back, cap/cover the GEM-containing tube strip and place on ice for no more than **1 h**.




Expose Wells at 45 Degrees



Transfer GEMs



GEMs

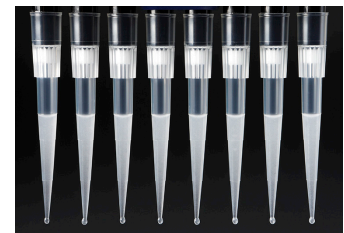

1.5 GEM-RT Incubation

Use a thermal cycler that can accommodate at least 100 μl volume. A volume of 125 μl is the preferred setting on the Bio-Rad C1000 Touch. In alternate thermal cyclers, use highest reaction volume setting.

- a. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
53°C	125 μl	~55 min

Step	Temperature	Time
1	53°C	00:45:00
2	85°C	00:05:00
3	4°C	Hold



- b. Store at 4°C for up to 72 h or at -20°C for up to a week, or proceed to the next step.

Step 2


Post GEM-RT Cleanup & cDNA Amplification

- 2.1 Post GEM-RT Cleanup – Dynabeads
- 2.2 cDNA Amplification
- 2.3 cDNA Cleanup – SPRIselect
- 2.4 cDNA QC & Quantification



2.0 Post GEM-RT Cleanup & cDNA Amplification



GET STARTED!				
Action	Item	10x PN	Preparation & Handling	Storage
Equilibrate to Room Temperature	Reducing Agent B	2000087	Thaw, vortex, verify no precipitate, centrifuge.	-20°C
	<input checked="" type="radio"/> Feature cDNA Primers 1 <i>Verify name & PN</i>	2000096	Vortex, centrifuge briefly.	-20°C
	 Beckman Coulter SPRIselect Reagent	-	Manufacturer's recommendations.	-
	Agilent Bioanalyzer High Sensitivity Kit If used for QC and quantification	-	Manufacturer's recommendations.	-
	Agilent TapeStation ScreenTape and Reagents If used for QC and quantification	-	Manufacturer's recommendations.	-
	Qubit dsDNA HS Assay Kit If used for QC and quantification	-	Manufacturer's recommendations.	-
	Dynabeads MyOne SILANE	2000048	Vortex thoroughly (≥30 sec) immediately before adding to the mix.	4°C
Place on ice	<input type="radio"/> Amp Mix	2000047/ 2000103	Vortex, centrifuge briefly.	-20°C
Thaw at 65°C	<input checked="" type="radio"/> Cleanup Buffer	2000088	Thaw for 10 min at 65°C at max speed on a thermomixer. Verify no visible crystals. Cool to room temperature.	-20°C
Obtain	<input type="radio"/> Recovery Agent	220016	-	Ambient
	Qiagen Buffer EB	-	Manufacturer's recommendations.	-
	Bio-Rad 10% Tween 20	-	Manufacturer's recommendations.	-
	10x Magnetic Separator	230003	-	Ambient
	Prepare 80% Ethanol Prepare 15 ml for 8 reactions.	-	-	-

2.1 Post GEM-RT Cleanup – Dynabeads

- a. Add **125 µl** Recovery Agent to each sample at room temperature. **DO NOT** pipette mix or vortex the biphasic mixture. Wait **60 sec**.

The resulting biphasic mixture contains Recovery Agent/Partitioning Oil (pink) and aqueous phase (clear), with no persisting emulsion (opaque).

If biphasic separation is incomplete, cap the tube strip, centrifuge briefly, and proceed to step b.




A smaller aqueous phase volume indicates a clog during GEM generation.



- b. Slowly remove **125 µl** Recovery Agent/Partitioning Oil (pink) from the bottom of the tube. **DO NOT** aspirate any aqueous sample.



- c. Prepare Dynabeads Cleanup Mix.

Dynabeads Cleanup Mix <i>Add reagents in the order listed</i>	PN	1X (µl)	4X + 10% (µl)	8X + 10% (µl)
● Cleanup Buffer	2000088	182	801	1602
 Dynabeads MyOne SILANE Vortex thoroughly (≥ 30 sec) immediately before adding to the mix. If still clumpy, pipette mix to resuspend completely. DO NOT centrifuge before use.	2000048	8	35	70
Reducing Agent B	2000087	5	22	44
Nuclease-free Water		5	22	44



- d. Vortex and add **200 µl** to each sample. Pipette mix 10x (pipette set to 200 µl).
- e. Incubate **10 min** at room temperature. Pipette mix again at **~5 min** after start of incubation to resuspend settled beads.



f. Prepare Elution Solution I. Vortex and centrifuge briefly.

Elution Solution I <i>Add reagents in the order listed</i>	PN	1X (μl)	10X (μl)
Buffer EB	-	98	980
10% Tween 20	-	1	10
<input type="radio"/> Reducing Agent B	2000084	1	10
Total	-	100	1000

g. At the end of **10 min** incubation, place on a 10x Magnetic Separator • **High** position (magnet • **High**) until the solution clears.

A white interface between the aqueous phase and Recovery Agent is normal.

h. Remove the supernatant.

i. Add **300 μl** 80% ethanol to the pellet while on the magnet. Wait **30 sec**.

j. Remove the ethanol.

k. Add **200 μl** 80% ethanol to pellet. Wait **30 sec**.

l. Remove the ethanol.

m. Centrifuge briefly. Place on the magnet • **Low**.n. Remove remaining ethanol. Air dry for **1 min**.o. Remove from the magnet. Immediately add **35.5 μl** Elution Solution I.

p. Pipette mix (pipette set to 30 μl) without introducing bubbles.

q. Incubate **2 min** at **room temperature**.r. Place on the magnet • **Low** until the solution clears.s. Transfer **35 μl** sample to a new tube strip.

2.2 cDNA Amplification



a. Prepare cDNA Amplification Mix on ice. Vortex and centrifuge briefly.

cDNA Amplification Reaction Mix <i>Add reagents in the order listed</i>	PN	1X (μl)	4X + 10% (μl)	8X + 10% (μl)
<input type="radio"/> Amp Mix	2000047/ 2000103	50	220	440
<input checked="" type="radio"/> Feature cDNA Primers 1 <i>Verify name & PN</i> <i>Use indicated primer only</i>	2000096	15	66	132
Total	-	65	286	572



b. Add **65 μl** cDNA Amplification Reaction Mix to **35 μl** sample.

c. Pipette mix 15x (pipette set to 90 μl). Centrifuge briefly.

d. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 μl	~30-45 min
Step	Temperature	Time
1	98°C	00:03:00
2	98°C	00:00:15
3	63°C <i>Version specific updated temperature</i>	00:00:20
4	72°C	00:01:00
5	Go to Step 2, see table below for total # of cycles	
6	72°C	00:01:00
7	4°C	Hold

Recommended starting point for cycle number optimization.


Cell Load	Total Cycles
<500	13
500-6,000	12
>6,000	11

The optimal number of cycles is a trade-off between generating sufficient final mass for library construction and minimizing PCR amplification artifacts. The number of cDNA cycles should also be reduced if large numbers of cells are sampled.




e. Store at **4°C** for up to **72 h** or proceed to the next step.


Step Overview
(steps 2.2 & 2.3)

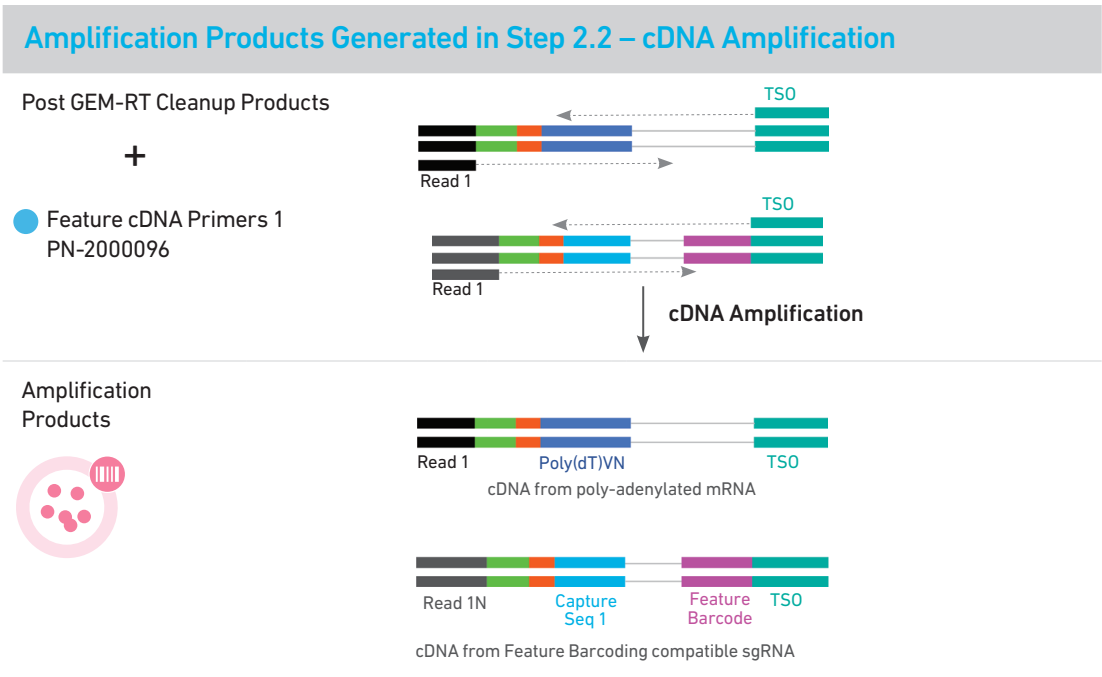

Step 2.3 – cDNA Cleanup – SPRIselect Overview

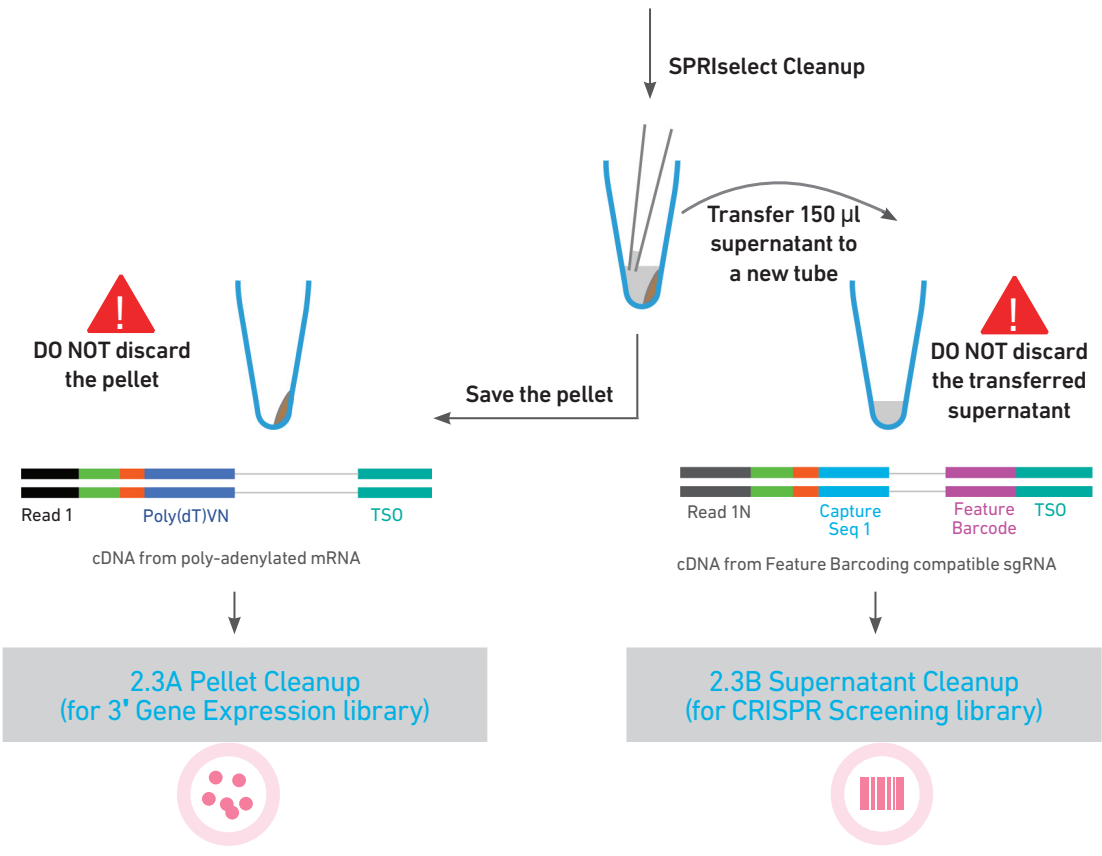


2.3 cDNA Cleanup – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add **60 µl** SPRIselect reagent (**0.6X**) to each sample and pipette mix 15x (pipette set to 150 µl).
- b. Incubate **5 min** at **room temperature**.
- c. Place on the magnet•**High** until the solution clears.
-  d. Transfer and save **150 µl** supernatant in a new tube strip without disturbing the pellet. Maintain at **room temperature**. **DO NOT** discard the transferred supernatant (cleanup for CRISPR Screening library construction).
- e. **DO NOT** discard the pellet (cleanup for 3' Gene Expression library construction). **Immediately** proceed to Pellet Cleanup (step 2.3A).



2.3A Pellet Cleanup


(for 3' Gene Expression library)

- i. Add **200 µl** 80% ethanol to the pellet. Wait **30 sec**.
- ii. Remove the ethanol.
- iii. **Repeat** steps i and ii for a total of 2 washes.
- iv. Centrifuge briefly and place on the magnet•**Low**.
- v. Remove any remaining ethanol. Air dry for **2 min**. **DO NOT** exceed **2 min** as this will decrease elution efficiency.
- vi. Remove from the magnet. Add **40.5 µl** Buffer EB. Pipette mix 15x.
- vii. Incubate **2 min** at **room temperature**.
- viii. Place the tube strip on the magnet•**High** until the solution clears.
- ix. Transfer **40 µl** sample to a new tube strip.
- x.  Store at **4°C** for up to **72 h** or at **-20°C** for up to **4 weeks**, or proceed to [step 2.4](#) followed by [step 3](#) for 3' Gene Expression Library Construction.



2.3B Transferred Supernatant Cleanup


(for CRISPR Screening library)

- i. Vortex to resuspend the SPRIselect reagent. Add **60 µl** SPRIselect reagent (**1.2X**) to **150 µl** of the transferred supernatant and pipette mix 15x (pipette set to 150 µl).
- ii. Incubate for **5 min** at **room temperature**.
- iii. Place on the magnet•**High** until the solution clears.
- iv. Remove supernatant.
- v. Add **200 µl** 80% ethanol to the pellet. Wait **30 sec**.
- vi. Remove the ethanol.
- vii. **Repeat** steps v and vi for a total of 2 washes.
- viii. Centrifuge briefly and place on the magnet•**Low**.
- ix. Remove any remaining ethanol. Air dry for **2 min**. **DO NOT** exceed **2 min** as this will decrease elution efficiency.
- x. Remove from the magnet. Add **50.5 µl** Buffer EB. Pipette mix 15x.
- xi. Incubate **2 min** at **room temperature**.
- xii. Place the tube strip on the magnet•**High** until the solution clears.
- xiii. Transfer **50 µl** sample to a new tube strip.
- xiv.  Store at **4°C** for up to **72 h** or at **-20°C** for up to **4 weeks**, or proceed directly to [step 4](#) for CRISPR Screening Library Construction.

2.4
Post cDNA Amplification
QC & Quantification

a. Run 1 μl of sample from **Pellet Cleanup** (step 2.3A-x), diluted 1:10 on an Agilent Bioanalyzer High Sensitivity chip. **DO NOT** run sample from 2.3B Transferred Supernatant Cleanup step.


For input cells with low RNA content (<1pg total RNA/cell), 1 μl undiluted product may be run. Lower molecular weight product (35 – 150 bp) may be present. This is normal and does not affect sequencing or application performance.

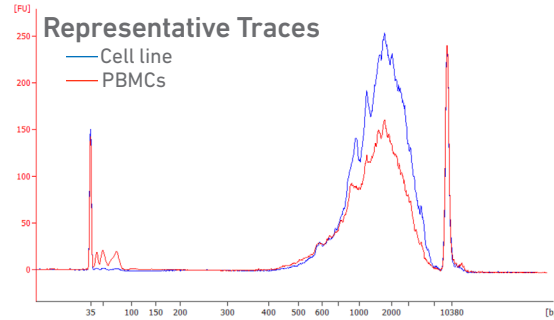


EXAMPLE CALCULATION

i. **Select Region**

Under the “Electropherogram” view choose the “Region Table”. Manually select the region of ~200 – ~9000 bp

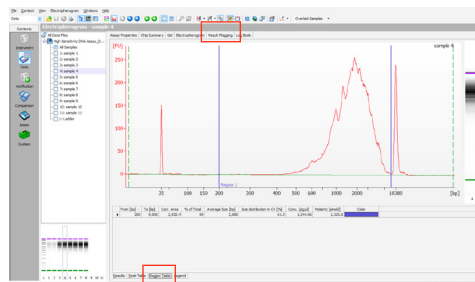

ii. **Note Concentration [pg/μl]**

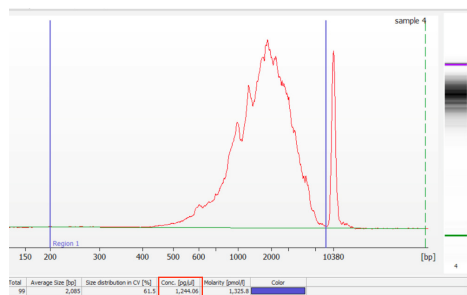

iii. **Calculate**

Multiply the cDNA concentration [pg/μl] reported via the Agilent 2100 Expert Software by the elution volume (40 μl) of the Post cDNA Amplification Reaction Clean Up sample (taking any dilution factors into account) and then divide by 1000 to obtain the total cDNA yield in ng.

Example Calculation of cDNA Total Yield

Concentration: 1244.06 pg/μl
Elution Volume: 40
Dilution Factor: 10

$$\begin{aligned} \text{Total cDNA Yield} &= \frac{\text{Conc'n (pg/}\mu\text{l}) \times \text{Elution Volume} \times \text{Dilution Factor}}{1000 \text{ (pg/ng)}} \\ &= \frac{1244.06(\text{pg}/\mu\text{l}) \times 40 \times 10}{1000 \text{ (pg/ng)}} \end{aligned}$$

! Carry forward **ONLY 25%** of total cDNA yield into 3' Gene Expression Library Construction (step 3)

VERSION SPECIFIC

$$\begin{aligned} &= 0.25 \times \text{Total cDNA yield} \\ &= 0.25 \times 497.62 = 124.4 \text{ ng} \end{aligned}$$

Refer to step 3.5 for appropriate number of Sample Index PCR cycles based on carry forward cDNA yield/input cDNA.

Alternate Quantification Methods:

- Agilent TapeStation. [See Appendix for representative traces](#)

Agilent Bioanalyzer or Agilent TapeStation are the recommended methods for accurate quantification.

(If using Qubit Fluorometer and Qubit dsDNA HS Assay Kit, [see Appendix](#))

Step 3

3' Gene Expression Library Construction

- 3.1 Fragmentation, End Repair & A-tailing
- 3.2 Post Fragmentation End Repair & A-tailing Double Sided Size Selection – SPRIselect
- 3.3 Adaptor Ligation
- 3.4 Post Ligation Cleanup – SPRIselect
- 3.5 Sample Index PCR
- 3.6 Post Sample Index PCR Double Sided Size Selection – SPRIselect
- 3.7 Post Library Construction QC

3.0 3' Gene Expression Library Construction



GET STARTED!				
Action	Item	10x PN	Preparation & Handling	Storage
Equilibrate to Room Temperature	● Fragmentation Buffer	2000091	Vortex, verify no precipitate, centrifuge briefly.	-20°C
	● Adaptor Oligos	2000094	Vortex, centrifuge briefly.	-20°C
	● Ligation Buffer	2000092	Vortex, verify no precipitate, centrifuge briefly.	-20°C
	● SI Primer	2000095	-	-20°C
	Chromium i7 Sample Index Plate	120262	-	-20°C
	Beckman Coulter SPRIselect Reagent	-	Manufacturer's recommendations.	-
	Agilent TapeStation Screen Tape and Reagents If used for QC		Manufacturer's recommendations.	
Agilent Bioanalyzer High Sensitivity kit If used for QC	-	Manufacturer's recommendations.	-	
Place on Ice	● Fragmentation Enzyme	2000090/ 2000104	Centrifuge briefly.	-20°C
	● DNA Ligase	220110/ 220131	Centrifuge briefly.	-20°C
	○ Amp Mix	2000047/ 2000103	Centrifuge briefly.	-20°C
	KAPA Library Quantification Kit for Illumina Platforms	-	Manufacturer's recommendations.	-
Obtain	Qiagen Buffer EB	-	-	Ambient
	10x Magnetic Separator	230003	See Tips & Best Practices.	Ambient
	Prepare 80% Ethanol Prepare 20 ml for 8 reactions	-	Prepare fresh.	Ambient

3.1 Fragmentation, End Repair & A-tailing

- a. Prepare a thermal cycler with the following incubation protocol.



Lid Temperature	Reaction Volume	Run Time
65°C	50 µl	~35 min
Step	Temperature	Time
Pre-cool block <i>Pre-cool block prior to preparing the Fragmentation Mix</i>	4°C	Hold
Fragmentation	32°C	00:05:00
End Repair & A-tailing	65°C	00:30:00
Hold	4°C	Hold

- b. Vortex Fragmentation Buffer. Verify there is no precipitate.

- c. Prepare Fragmentation Mix on ice. Pipette mix and centrifuge briefly.

Fragmentation Mix <i>Add reagents in the order listed</i>	PN	1X (µl)	4X + 10% (µl)	8X + 10% (µl)
● Fragmentation Buffer	2000091/ 2000104	5	22	44
● Fragmentation Enzyme	2000090	10	44	88
Total	-	15	66	132



- d. Transfer **ONLY 10 µl** purified cDNA sample from Pellet Cleanup (step 2.3A-x) to a tube strip.

Note that only **10 µl** (25%) cDNA sample transfer is sufficient for generating 3' Gene Expression library.

The remaining **30 µl** (75%) cDNA sample can be stored at **4°C** for up to **72 h** or at **-20°C** for up to **4 weeks** for generating additional 3' Gene Expression libraries.

- e. Add **25 µl** Buffer EB to each sample.

- f. Add **15 µl** Fragmentation Mix to each sample.

- g. Pipette mix 15x (pipette set to 35 µl) on ice. Centrifuge briefly.

- h. Transfer into the pre-cooled thermal cycler (**4°C**) and press "SKIP" to initiate the protocol.

3.2
Post Fragmentation,
End Repair & A-tailing
Double Sided Size
Selection – SPRIselect

- a. Vortex to resuspend SPRIselect reagent. Add **30 μ l** SPRIselect (**0.6X**) reagent to each sample. Pipette mix 15x (pipette set to 75 μ l).
- b. Incubate **5 min** at **room temperature**.
- c. Place on the magnet•**High** until the solution clears. **DO NOT** discard supernatant.



- d. Transfer **75 μ l** supernatant to a new tube strip.
- e. Vortex to resuspend SPRIselect reagent. Add **10 μ l** SPRIselect reagent (**0.8X**) to each sample. Pipette mix 15x (pipette set to 80 μ l).
- f. Incubate **5 min** at **room temperature**.
- g. Place on the magnet•**High** until the solution clears.



- h. Remove **80 μ l** supernatant. **DO NOT** discard any beads.
- i. Add **125 μ l** 80% ethanol to the pellet. Wait **30 sec**.
- j. Remove the ethanol.
- k. **Repeat** steps i and j for a total of 2 washes.
- l. Centrifuge briefly. Place on the magnet•**Low** until the solution clears. Remove remaining ethanol. **DO NOT** over dry to ensure maximum elution efficiency.
- m. Remove from the magnet. Add **50.5 μ l** Buffer EB to each sample. Pipette mix 15x.
- n. Incubate **2 min** at **room temperature**.
- o. Place on the magnet•**High** until the solution clears.
- p. Transfer **50 μ l** sample to a new tube strip.

3.3 Adaptor Ligation

a. Prepare Adaptor Ligation Mix. Pipette mix and centrifuge briefly.

Adaptor Ligation Mix <i>Add reagents in the order listed</i>	PN	1X (µl)	4X + 10% (µl)	8X + 10% (µl)
● Ligation Buffer	2000092	20	88	176
● DNA Ligase	220110/ 220131	10	44	88
● Adaptor Oligos	2000094	20	88	176
Total	-	50	220	440

b. Add 50 µl Adaptor Ligation Mix to 50 µl sample. Pipette mix 15x (pipette set to 90 µl). Centrifuge briefly.

c. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
30°C	100 µl	15 min
Step	Temperature	Time
1	20°C	00:15:00
2	4°C	Hold

3.4 Post Ligation Cleanup – SPRIselect

- a. Vortex to resuspend SPRIselect Reagent. Add **80 µl** SPRIselect Reagent (**0.8X**) to each sample. Pipette mix 15x (pipette set to 150 µl).
- b. Incubate **5 min** at **room temperature**.
- c. Place on the magnet•**High** until the solution clears.
- d. Remove the supernatant.
- e. Add **200 µl** 80% ethanol to the pellet. Wait **30 sec**.
- f. Remove the ethanol.
- g. **Repeat** steps e and f for a total of 2 washes.
- h. Centrifuge briefly. Place on the magnet•**Low**.
- i. Remove any remaining ethanol. Air dry for **2 min**.
DO NOT exceed **2 min** as this will decrease elution efficiency.
- j. Remove from the magnet. Add **30.5 µl** Buffer EB. Pipette mix 15x.
- k. Incubate **2 min** at **room temperature**.
- l. Place on the magnet•**Low** until the solution clears.
- m. Transfer **30 µl** sample to a new tube strip.

3.5 Sample Index PCR



a. Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run. Record the 10x Sample Index name (PN-220103 Chromium i7 Sample Index Plate well ID) used.

b. Prepare Sample Index PCR Mix.

Sample Index PCR Mix <i>Add reagents in the order listed</i>	PN	1X (μl)	4X + 10% (μl)	8X + 10% (μl)
<input type="radio"/> AmpMix	2000047/ 2000103	50	220	440
<input checked="" type="radio"/> SI Primer	2000095	10	44	88
Total	-	60	264	528

c. Add 60 μl Sample Index PCR Mix to 30 μl sample.

d. Add 10 μl of an individual Chromium i7 Sample Index to each well and record the well ID used. Pipette mix 5x (pipette set to 90 μl). Centrifuge briefly.

e. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 μl	~25-40 min
Step	Temperature	Time
1	98°C	00:00:45
2	98°C	00:00:20
3	54°C	00:00:30
4	72°C	00:00:20
5	Go to step 2, see below for # of cycles	
6	72°C	00:01:00
7	4°C	Hold



The total cycles should be optimized based on 25% carry forward cDNA yield/input calculated during cDNA QC & Quantification (step 2.4)

Recommended cycle numbers

cDNA Input	Total Cycles
1-25 ng	14-16
25-150 ng	12-14
150-500 ng	10-12
500-1,000 ng	8-10
1,000-1,500 ng	6-8



f. Store at 4°C for up to 72 h or proceed to the next step.


3.6
Post Sample Index
PCR Double Sided Size
Selection – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add **60 µl** SPRIselect Reagent (**0.6X**) to each sample. Pipette mix 15x (pipette set to 150 µl).
- b. Incubate **5 min** at **room temperature**.
- c. Place the magnet•**High** until the solution clears. **DO NOT** discard supernatant.
- d. Transfer **150 µl** supernatant to a new tube strip.
- e. Vortex to resuspend the SPRIselect reagent. Add **20 µl** SPRIselect Reagent (**0.8X**) to each sample. Pipette mix 15x (pipette set to 150 µl).
- f. Incubate **5 min** at **room temperature**.
- g. Place the magnet•**High** until the solution clears.
- h. Remove **165 µl** supernatant. **DO NOT** discard any beads.
- i. With the tube still in the magnet, add **200 µl** 80% ethanol to the pellet. Wait **30 sec**.
- j. Remove the ethanol.
- k. **Repeat** steps i and j for a total of 2 washes.
- l. Centrifuge briefly. Place on the magnet•**Low**. Remove remaining ethanol.
- m. Remove from the magnet. Add **35.5 µl** Buffer EB. Pipette mix 15x.
- n. Incubate **2 min** at **room temperature**.
- o. Place on the magnet•**Low** until the solution clears.
- p. Transfer **35 µl** to a new tube strip.
- q. Store at **4°C** for up to **72 h** or at **-20°C** for **long-term** storage.



3.7 Post Library Construction QC

Run 1 μl sample at 1:10 dilution on an Agilent Bioanalyzer High Sensitivity chip.

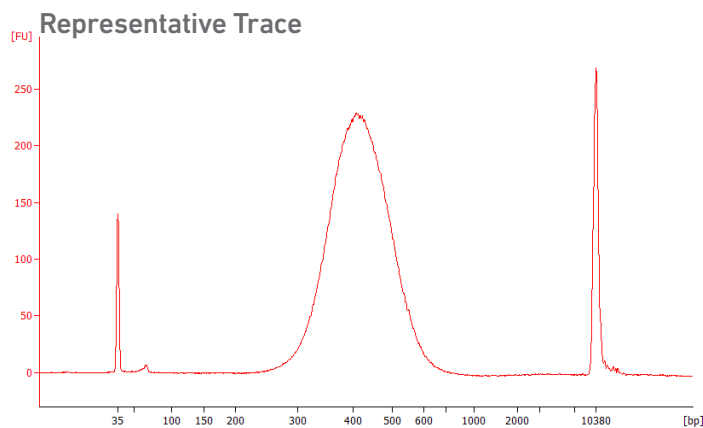

Determine the average fragment size from the Bioanalyzer trace. This will be used as the insert size for library quantification.

Alternate QC Method:

- Agilent TapeStation. [See Appendix for representative traces](#)

[See Appendix for Post Library Construction Quantification](#)

Step 4

CRISPR Screening Library Construction

- 4.1 Guide RNA cDNA Cleanup– SPRIselect
- 4.2 Feature PCR
- 4.3 Post Feature PCR Cleanup – SPRIselect
- 4.4 Sample Index PCR
- 4.5 Post Sample Index PCR Double Sided Size Selection – SPRIselect
- 4.6 Post Library Construction QC

4.0 CRISPR Screening Library Construction



GET STARTED!				
Action	Item	10x PN	Preparation & Handling	Storage
Equilibrate to Room Temperature	● Feature SI Primers 1	2000098	-	-20°C
	● Feature SI Primers 2	2000099	-	-20°C
	Chromium i7 Sample Index Plate	120262	-	-20°C
	Beckman Coulter SPRIselect Reagent	-	Manufacturer's recommendations.	-
	Agilent TapeStation Screen Tape and Reagents If used for QC	-	Manufacturer's recommendations.	-
	Agilent Bioanalyzer High Sensitivity kit If used for QC	-	Manufacturer's recommendations.	-
Place on Ice	○ Amp Mix	2000047	Centrifuge briefly.	-20°C
	KAPA Library Quantification Kit for Illumina Platforms	-	Manufacturer's recommendations.	-
Obtain	Qiagen Buffer EB	-	-	Ambient
	10x Magnetic Separator	230003	See Tips & Best Practices.	Ambient
	Prepare 80% Ethanol Prepare 20 ml for 8 reactions	-	Prepare fresh.	Ambient

4.1 Guide RNA cDNA Cleanup – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add **50 µl** SPRIselect reagent (**1.0X**) to **50 µl** Transferred Supernatant Cleanup (step 2.3B-xiv) and pipette mix 15x (pipette set to 150 µl).
- b. Incubate **5 min** at **room temperature**.
- c. Place on the magnet•**High** until the solution clears.
- d. Remove supernatant.
- e. Add **200 µl** 80% ethanol to the supernatant. Wait **30 sec**.
- f. Remove the ethanol.
- g. **Repeat** steps e and f for a total of 2 washes.
- h. Centrifuge briefly and place on the magnet•**Low**.
- i. Remove any remaining ethanol. Air dry for **2 min**. **DO NOT** exceed **2 min** as this will decrease elution efficiency.
- j. Remove from the magnet. Add **30.5 µl** Buffer EB. Pipette mix 15x.
- k. Incubate **2 min** at **room temperature**.
- l. Place the tube strip on the magnet•**High** until the solution clears.
- m. Transfer **30 µl** sample to a new tube strip.
- n. Store at **4°C** for up to **72 h** or at **-20°C** for up to a **week**, or proceed to the next step.



4.2 Feature PCR

a. Prepare Feature PCR Mix on ice. Vortex and centrifuge briefly.

Feature PCR Mix <i>Add reagents in the order listed</i>	PN	1X (μl)	4X + 10% (μl)	8X + 10% (μl)
○ Amp Mix	2000047	50	220	440
● Feature SI Primers 1	2000098	45	198	396
Total	-	95	418	836

b. Transfer **ONLY 5 μl** from Guide RNA cDNA Cleanup (step 4.1m) to a new tube strip.

Note that only **5 μl** of the DNA sample transfer is sufficient for generating CRISPR Screening library. The remaining **25 μl** sample can be stored at **4°C** for up to **72 h** or at **-20°C** for up to **4 weeks**, for generating additional CRISPR Screening libraries.

c. Add **95 μl** Feature PCR Mix to **5 μl** sample.

d. Pipette mix 15x (pipette set to 90 μl). Centrifuge briefly.

e. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 μl	~30-45 min
Step	Temperature	Time
1	98°C	00:00:45
2	98°C	00:00:20
3	60°C	00:00:05
4	72°C	00:00:05
5	Go to Step 2, repeat 14X for a total of 15 cycles	
6	72°C	00:01:00
7	4°C	Hold

4.3 Post Feature PCR Cleanup – SPRIselect

- a. Vortex to resuspend SPRIselect Reagent. Add **100 μ l** SPRIselect Reagent (**1.0X**) to each sample. Pipette mix 15x (pipette set to 150 μ l).
- b. Incubate **5 min** at **room temperature**.
- c. Place on the magnet•**High** until the solution clears.
- d. Remove the supernatant.
- e. Add **300 μ l** 80% ethanol to the pellet. Wait **30 sec**.
- f. Remove the ethanol.
- g. **Repeat** steps e and f for a total of 2 washes.
- h. Centrifuge briefly. Place on the magnet•**Low**.
- i. Remove any remaining ethanol. Air dry for **1 min**.
DO NOT exceed **1 min** as this will decrease elution efficiency.
- j. Remove from the magnet. Add **30.5 μ l** Buffer EB. Pipette mix 15x.
- k. Incubate **2 min** at **room temperature**.
- l. Place on the magnet•**Low** until the solution clears.
- m. Transfer **30 μ l** sample to a new tube strip.

4.4 Sample Index PCR



a. Choose the appropriate sample index sets to ensure that no sample indices overlap in a multiplexed sequencing run. Avoid using the same sample index sets for both 3' Gene Expression and CRISPR Screening libraries constructed from the same sample. Record the 10x Sample Index name (PN-220103 Chromium i7 Sample Index Plate well ID) used.

b. Prepare Sample Index PCR Mix.

Sample Index PCR Mix <i>Add reagents in the order listed</i>	PN	1X (μl)	4X + 10% (μl)	8X + 10% (μl)
<input type="radio"/> AmpMix	2000047	50	220	440
<input checked="" type="radio"/> Feature SI Primers 2	2000099	35	154	308
Total	-	85	374	748

c. Transfer **ONLY 5 μl** Post Feature PCR Cleanup sample (step 4.3m) to a new tube strip. Note that only **5 μl** sample transfer is sufficient for generating CRISPR Screening library. The remaining **25 μl** sample can be stored at **4°C** for up to **72 h** or at **-20°C** for up to **4 weeks**, for generating additional CRISPR Screening libraries.

d. Add **85 μl** Sample Index PCR Mix to **5 μl** sample (Post Feature PCR Cleanup).

e. Add **10 μl** of an individual Chromium i7 Sample Index to each well and record their assignment. Pipette mix 5x (pipette set to 90 μl). Centrifuge briefly.

f. Incubate in a thermal cycler with the following protocol.

Lid Temperature	Reaction Volume	Run Time
105°C	100 μl	~25-40 min
Step	Temperature	Time
1	98°C	00:00:45
2	98°C	00:00:20
3	54°C	00:00:30
4	72°C	00:00:20
5	Go to step 2, repeat 4X for a total of 5 cycles	
6	72°C	00:01:00
7	4°C	Hold


4.5
Post Sample Index
PCR Double Sided Size
Selection – SPRIselect

- a. Vortex to resuspend the SPRIselect reagent. Add **70 μ l** SPRIselect Reagent (**0.7X**) to each sample. Pipette mix 15x (pipette set to 150 μ l).
- b. Incubate **5 min** at **room temperature**.
- c. Place the magnet•**High** until the solution clears. **DO NOT** discard supernatant.
- d. Transfer **160 μ l** supernatant to a new tube strip.
- e. Vortex to resuspend the SPRIselect reagent. Add **30 μ l** SPRIselect Reagent (**1.0X**) to each sample. Pipette mix 15x (pipette set to 150 μ l).
- f. Incubate **5 min** at **room temperature**.
- g. Place the magnet•**High** until the solution clears.
- h. Remove the supernatant.
- i. Add **300 μ l** 80% ethanol to the pellet. Wait **30 sec**.
- j. Remove the ethanol.
- k. Add **200 μ l** 80% ethanol to the pellet. Wait **30 sec**.
- l. Remove the ethanol.
- m. Centrifuge briefly. Place on the magnet•**Low**.
- n. Remove remaining ethanol. Air dry for 1 min.
- o. Remove from the magnet. Add **30.5 μ l** Buffer EB. Pipette mix 15x.
- p. Incubate **2 min** at **room temperature**.
- q. Place on the magnet•**Low** until the solution clears.
- r. Transfer **30 μ l** to a new tube strip.
- s. Store at **4°C** for up to **72 h** or at **-20°C** for **long-term** storage.



4.6 Post Library Construction QC

Run 1 μl sample at 1:25 dilution on an Agilent Bioanalyzer High Sensitivity chip.

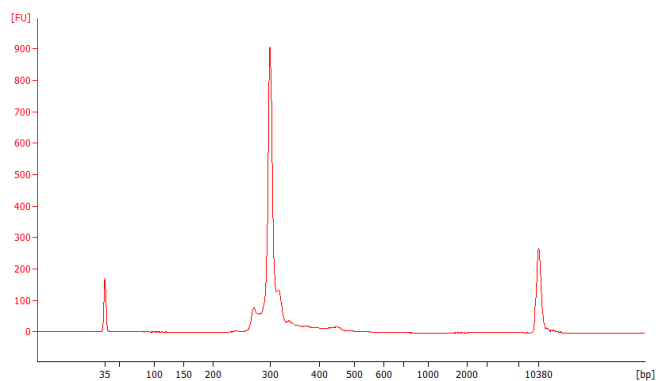

Determine the average fragment size from the Bioanalyzer trace. This will be used as the insert size for library quantification.

Alternate QC Method:

- Agilent TapeStation. [See Appendix for representative traces](#)

[See Appendix for Post Library Construction Quantification](#)


Sequencing

5


Sequencing Libraries

Chromium Single Cell 3' Gene Expression and CRISPR Screening libraries comprise standard Illumina paired-end constructs which begin with P5 and end with P7. These libraries include 16 bp 10x Barcodes at the start of TruSeq Read 1 and Nextera Read 1 (Read 1N) respectively while 8 bp sample index sequences are incorporated as the i7 index read. TruSeq Read 1 and TruSeq Read 2 are standard Illumina sequencing primer sites used in paired-end sequencing of Single Cell 3' Gene Expression libraries. Nextera Read 1 (Read 1N) and TruSeq Read 2 are used for paired-end sequencing of Single Cell 3' CRISPR Screening libraries. Sequencing these libraries produce a standard Illumina BCL data output folder.

Chromium Single Cell 3' Gene Expression Library

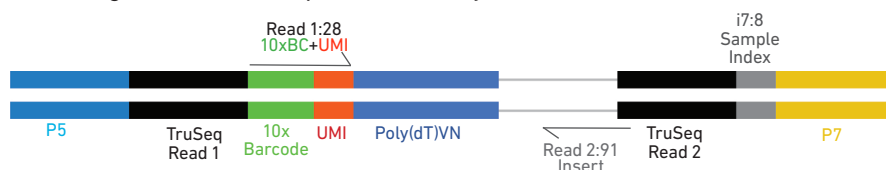

Chromium Single Cell 3' CRISPR Screening Library

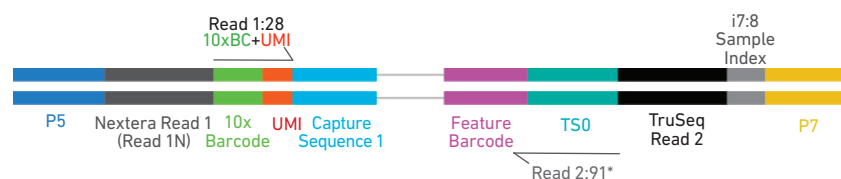

*Minimum required Read 2 length for CRISPR Screening libraries is 70 bp

Illumina Sequencer Compatibility

The compatibility of the listed sequencers has been verified by 10x Genomics. Some variation in assay performance is expected based on sequencer choice. For more information about performance variation, visit the 10x Genomics Support website.

- MiSeq
- NextSeq 500/550*
- HiSeq 2500 (Rapid Run)
- HiSeq 3000/4000
- NovaSeq

*Sequencing Chromium Single Cell libraries on the NextSeq 500/550 platform may yield reduced sequence quality and sensitivity relative to the MiSeq, HiSeq, and NovaSeq platforms. Refer to the 10x Genomics Support website for more information.

Sample Indices

Each sample index in the Chromium i7 Sample Index Kit (PN-120262) is a mix of 4 different sequences to balance across all 4 nucleotides. If multiple samples are pooled in a sequence lane, the sample index name (i.e. the Chromium i7 Sample Index plate well ID, SI-GA-) is needed in the sample sheet used for generating FASTQs with "cellranger mkfastq". Samples utilizing the same sample index should not be pooled together, or run on the same flow cell lane, as this would not enable correct sample demultiplexing.

3' Gene Expression Library Sequencing Depth & Run Parameters

Sequencing Depth	Minimum 20,000 read pairs per cell
Sequencing Type	Paired-end, single indexing
Sequencing Read	Recommended Number of Cycles
Read 1	28 cycles
i7 Index	8 cycles
i5 Index	0 cycles
Read 2	91 cycles

CRISPR Screening Library Sequencing Depth & Run Parameters

Sequencing Depth	Minimum 5,000 read pairs per cell
Sequencing Type	Paired-end, single indexing
Sequencing Read	Recommended Number of Cycles
Read 1	28 cycles
i7 Index	8 cycles
i5 Index	0 cycles
Read 2	91 cycles

Library Loading

Once quantified and normalized, the 3' Gene Expression and CRISPR Screening libraries should be denatured and diluted as recommended for Illumina sequencing platforms. Refer to Illumina documentation for denaturing and diluting libraries. Refer to the 10x Genomics Support website, for more information.

Instrument	3' Gene Expression libraries only or 3' Gene Expression + CRISPR Screening libraries		CRISPR Screening libraries only	
	Loading Concentration (pM)	PhiX (%)	Loading Concentration (pM)	PhiX (%)
MiSeq	11	1	11	1
NextSeq 500/550	1.8	1	1.8	20**
HiSeq 2500 (RR)	11	1	11	1
HiSeq 4000	240	1	240	1
NovaSeq	150*/300	1	150*/300	1

* Use 150pM loading concentration for Illumina XP workflow.

** When sequencing CRISPR screening libraries alone to create more diverse clusters for 30 cycles of Read 2.

Library Pooling

The 3' Gene Expression and CRISPR Screening libraries maybe pooled for sequencing, taking into account the differences in cell number and per-cell read depth requirements between each library. Samples utilizing the same sample index should not be pooled together, or run on the same flow cell lane, as this would not enable correct sample demultiplexing.

Library Pooling Example:

Libraries	Sequencing Depth (read pairs per cell)	Library Pooling Ratio
3' Gene Expression library	20,000	4
CRISPR Screening library	5,000	1

Troubleshooting



6.1 GEMs

STEP NORMAL REAGENT CLOGS & WETTING FAILURES

1.4 d
After Chip B is removed from the Controller and the wells are exposed



All 8 Recovery Wells are similar in volume and opacity.




Recovery Well G indicates a reagent clog. Recovery Well C and E indicate a wetting failure. Recovery Wells B, D, and F are normal. Wells A and H contain 50% Glycerol Solution.


The image indicates clogs in the Gel Bead line (orange arrow) and the sample line (yellow arrow) as evidenced by higher than usual volumes in the input wells.



1.4 f
Transfer GEMs from Chip E Recovery Wells

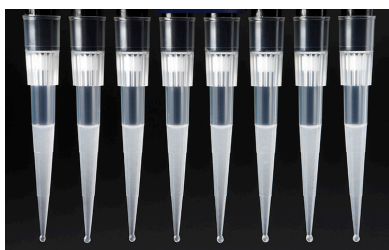


All liquid levels are similar in volume and opacity without air trapped in the pipette tips.

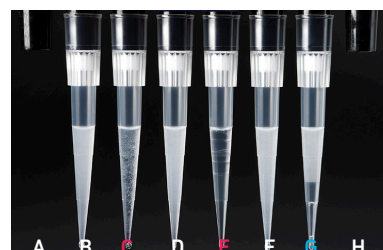


Pipette tips C and E indicate a wetting failure. Pipette tip C contains partially emulsified GEMs. Emulsion is absent in pipette tip E. Pipette tip G indicates a reagent clog.

STEP **NORMAL** **REAGENT CLOGS & WETTING FAILURES**

2.1 a
After transfer of the
GEMs +
Recovery Agent

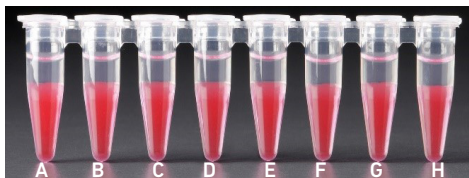

All liquid levels are similar in the aqueous sample volume (clear) and Recovery Agent/Partitioning Oil (pink).

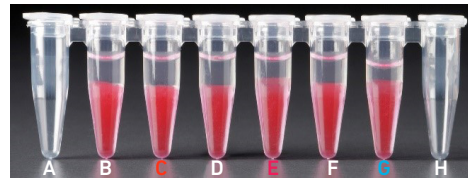

Tube **G** indicates a reagent clog has occurred. There is a decreased volume of aqueous layer (clear).
Tube **C** and **E** indicate a wetting failure has occurred. There is an abnormal residual volume of Recovery Agent/Partitioning Oil (pink).

2.1 b
After aspiration of
Recovery Agent/
Partitioning Oil



All liquid volumes are similar in the aqueous sample volume (clear) and residual Recovery Agent/Partitioning Oil (pink).



Tube **G** indicates a reagent clog has occurred. There is a decreased volume of aqueous layer (clear). There is also a greater residual volume of Recovery Agent/Partitioning Oil (pink).
Tube **C** and **E** indicate a wetting failure has occurred. There is an abnormal residual volume of Recovery Agent/Partitioning Oil (pink).

2.1 d
After addition of
Dynabeads
Cleanup Mix



All liquid volumes are similar after addition of the Dynabeads Cleanup Mix.



Tube **G** indicates a reagent clog has occurred. There is an abnormal ratio of Dynabeads Cleanup Mix (brown) to Recovery Agent/Partitioning Oil (appears white).
Tube **C** and **E** indicate a wetting failure has occurred. There is an abnormal ratio of Dynabeads Cleanup Mix (brown) to Recovery Agent/Partitioning Oil (appears white).

If a channel clogs or wetting failure occurs during GEM generation, it is recommended that the sample be remade. If any of the listed issues occur, take a picture and send it to support@10xgenomics.com for further assistance.

6.2 Chromium Controller Errors

If the Chromium Controller or the Chromium Single Cell Controller fails to start, an error tone will sound and one of the following error messages will be displayed:

- a. **Chip not read – Try again:** Eject the tray, remove and/or reposition the 10x Chip Holder assembly and try again. If the error message is still received after trying this more than twice, contact support@10xgenomics.com for further assistance.
- b. **Check gasket:** Eject the tray by pressing the eject button to check there is a 10x Gasket on the Chromium Chip. In the case when the 10x Gasket installation was forgotten, install and try again. In the case when a 10x Gasket was already installed, remove, reapply, and try again. If the error message is still received after trying either of these more than twice, contact support@10xgenomics.com for further assistance.
- c. **Pressure not at Setpoint:**
 - i. If this message is received within a few seconds of starting a run, eject the tray by pressing the eject button and check for dirt or deposits on the 10x Gasket. If dirt is observed, replace with a new 10x Gasket and try again. If the error message is still received after trying this more than twice, contact support@10xgenomics.com for further assistance.
 - ii. If this message is received after a few minutes into the run, the Chromium Chip must be discarded. **Do not try running this Chromium Chip again as this may damage the Chromium Controller.**
- d. **CAUTION: Chip Holder not Present:** Eject the tray by pressing the eject button to check there is a 10x Chip Holder encasing the Chromium Chip. In the case when the 10x Chip Holder was forgotten, install with a 10x Gasket in place, and try again. If the error message is still received after a 10x Chip Holder is confirmed as in place, contact support@10xgenomics.com for further assistance.
- e. **Invalid Chip CRC Value:** This indicates the Chromium Chip has encountered an error, should not be run, and must be discarded. Contact support@10xgenomics.com for further assistance.

Appendix

Post Library Construction Quantification

Agilent TapeStation Traces

Compatible sgRNA Specifications

Oligonucleotide Sequences

7

Post Library Construction Quantification

- a. Thaw KAPA Library Quantification Kit for Illumina Platforms.
- b. Dilute **2 µl** sample with deionized water to appropriate dilutions that fall within the linear detection range of the KAPA Library Quantification Kit for Illumina Platforms. (For more accurate quantification, make the dilution(s) in duplicate).
- c. Make enough Quantification Master Mix for the DNA dilutions per sample and the DNA Standards (plus 10% excess) using the guidance for 1 reaction volume below.

Quantification Master Mix	1X (µl)
SYBR Fast Master Mix + Primer	12
Water	4
Total	16

- d. Dispense **16 µl** Quantification Master Mix for sample dilutions and DNA Standards into a 96 well PCR plate.
- e. Add **4 µl** sample dilutions and **4 µl** DNA Standards to appropriate wells. Centrifuge briefly.
- f. Incubate in a thermal cycler with the following protocol.

Step	Temperature	Run Time
1	95°C	00:03:00
2	95°C	00:00:05
3	67°C	00:00:30
4	Go to Step 2, 29X (Total 30 cycles)	


- g. Follow the manufacturer's recommendations for qPCR-based quantification. For library quantification for sequencer clustering, determine the concentration based on insert size derived from the Bioanalyzer/TapeStation trace.

Agilent TapeStation Traces


Agilent TapeStation Traces

Agilent TapeStation High Sensitivity D5000 ScreenTape was used. Protocol steps correspond to the Chromium Single Cell 3 v3 Reagent Kits User Guide with Feature Barcoding technology for CRISPR Screening (CG000184)

Protocol Step 2.4 – cDNA QC & Quantification

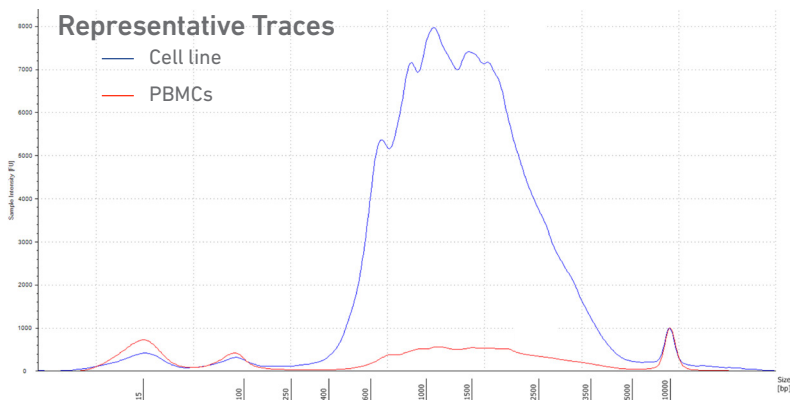

Protocol Step 4.6 – Post Library Construction QC

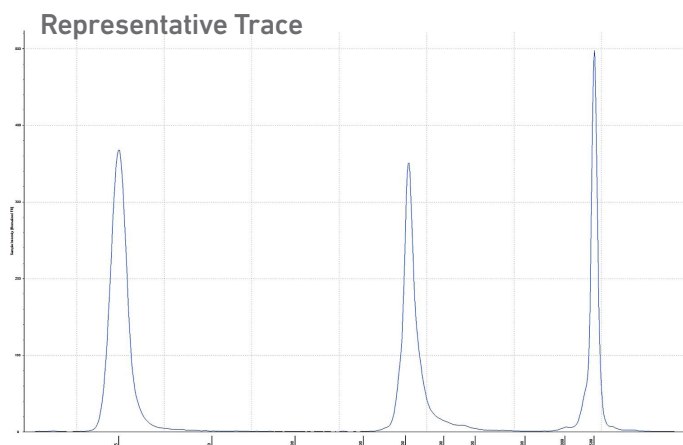

Alternate QC Method:

[Qubit Fluorometer and Qubit dsDNA HS Assay Kit](#)

Multiply the cDNA concentration reported via the Qubit Fluorometer by the elution volume (40 μ l) to obtain the total cDNA yield in ng. To determine the equivalent range using the Agilent 2100 Expert Software, select the region encompassing 35-10,000 bp.

Compatible sgRNA Specifications


Integration of Capture Sequence 1 and Capture Sequence 2 in sgRNA

A representative sgRNA sequence along with the specific capture sequences integrated in two different locations in the sgRNA are shown

Capture Sequence 1


Capture Sequence 1 on Gel Bead: 5'-TTGCTAGGACCGCCTTAAAGC-3'

Capture Sequence 1 integrated in sgRNA hairpin

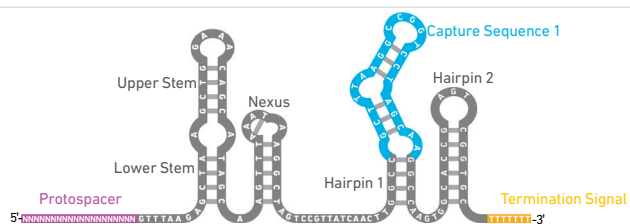

5'-NNNNNNNNNNNNNNNNNNNNNGTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTggccGCTTTAAGGCCGGTCTAGCAAggccAAGTGGCACCGAGTCGGTGC TTTTTT-3'

Capture Sequence 1 integrated in sgRNA 3'-end

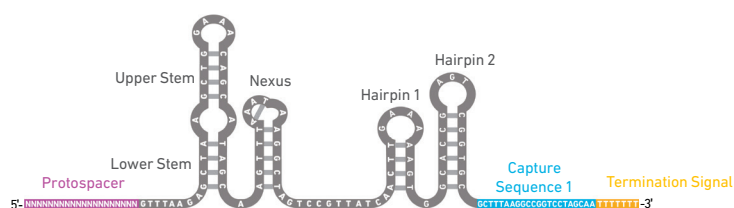


5'-NNNNNNNNNNNNNNNNNNNNNGTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTgaaaAAGTGGCACCGAGTCGGTGC GCTTTAAGGCCGGTCTAGCAA TTTTTT-3'

Capture Sequence 2


Capture Sequence 2 on Gel Bead oligonucleotide strand: 5'-CCTTAGCCGCTAATAGGTGAGC-3'

Capture Sequence 2 integrated in sgRNA hairpin

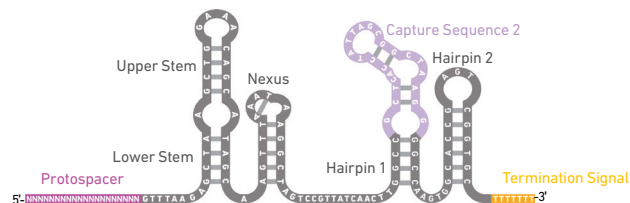

5'-NNNNNNNNNNNNNNNNNNNNNGTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTggccGCTCACCTATTAGCGGCTAAGGggccAAGTGGCACCGAGTCGGTGC TTTTTT-3'

Capture Sequence 2 integrated in sgRNA 3'-end

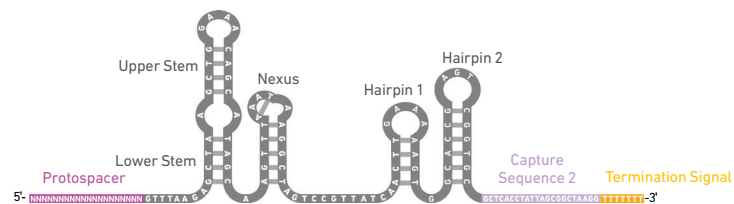

5'-NNNNNNNNNNNNNNNNNNNNNGTTTAAGAGCTAAGCTGGAAACAGCATAGCAAGTTTAAATAAGGCTAGTCCGTTATCAACTTgaaaAAGTGGCACCGAGTCGGTGC GCTCACCTATTAGCGGCTAAGG TTTTTT-3'


Oligonucleotide Sequences

Protocol steps correspond to the Chromium Single Cell 3 v3 Reagent Kits User Guide with Feature Barcoding technology for CRISPR Screening (CG000184)

Protocol Step 1.5 – GEM-RT Incubation

Gel Bead
Primers

5'-CTACACGACGCTCTCCGATCT-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-TTTTTTTTTTTTTTTTTTTTTTTTTTTT-3'

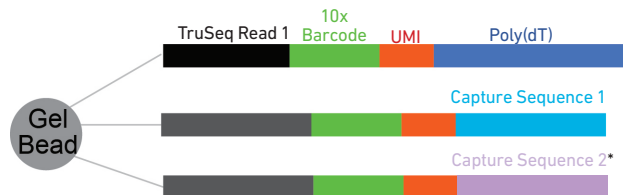

*Capture Sequence 2
is not illustrated in the
protocol steps below

5'-GTCAGATGTGTATAAGAGACAG-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-TTGCTAGGACCGCCCTAAAGC-3'

Template Switch
Oligonucleotide



TSO

5'-AAGCAGTGGTATCAACGCAGAGTACATrGrGrG-3'

PCR Products

cDNA from poly-adenylated mRNA



5'-CTACACGACGCTCTCCGATCT-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-TTTTTTTTTTTTTTTTTTTTTTTTTTTT-cDNA_Insert-CCCATGTACTCTGCGTTGATACCACTGCTT-3'

cDNA from sgRNA



5'-GTCAGATGTGTATAAGAGACAG-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-TTGCTAGGACCGCCCTAAAGC-sgRNA_Feature_Barcode-CCCATGTACTCTGCGTTGATACCACTGCTT-3'

Protocol Step 2.2 – cDNA Amplification

Feature cDNA
Primers 1
PN-2000096

[Amplifies cDNA from poly-adenylated mRNA](#)

Forward Primer: Partial Read 1

5'-CTACACGACGCTCTCCGATCT-3'

Reverse Primer: Partial TSO

5'-AAGCAGTGGTATCAACGCAGAG-3'

[Amplifies cDNA from sgRNA](#)

Forward Primer: Partial Read 1N


5'-GCAGCGTCAGATGTGTATAAGAGACAG-3'

Reverse primer: Partial TSO

5'-AAGCAGTGGTATCAACGCAGAG-3'


Amplification
Products

Amplified cDNA from poly-adenylated mRNA

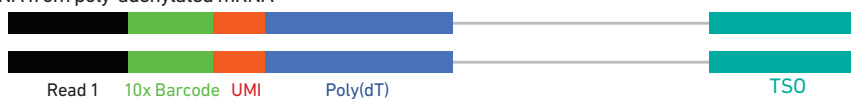

5'-CTACACGACGCTCTCCGATCT-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-TTTTTTTTTTTTTTTTTTTTTTTTTTTT-cDNA_Insert-CCCATGTACTCTGCGTTGATACCACTGCTT-3'
3'-GATGTCTGCGAGAAGGCTAGA-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-AAAAAAAAAAAAAAAAAAAAAAAAAAAA-cDNA_Insert-GGGTACATGAGACGCAACTATGGTGACGAA-5'


Amplified cDNA from sgRNA

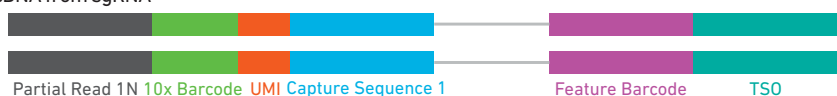


5'-GCAGCGTCAGATGTGTATAAGAGACAG-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-TTGCTAGGACCGCCCTAAAGC-sgRNA_Feature_Barcode-CCCATGTACTCTGCGTTGATACCACTGCTT-3'
3'-CGTCGAGTCTACACATATTCTGTCT-NNNNNNNNNNNNNNNN-NNNNNNNNNNNN-AACGATCTGGCCGGAATTTCG-sgRNA_Feature_Barcode-GGGTACATGAGACGCAACTATGGTGACGAA-5'

Protocol Step 4.2 – Feature PCR (for CRISPR Screening Library Construction)

Feature SI
Primers 2
PN-2000099

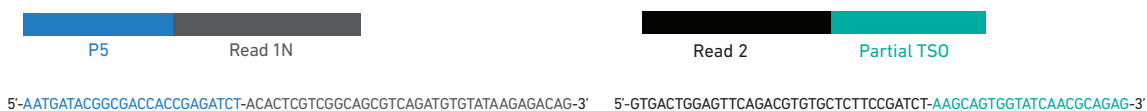


Feature PCR Product

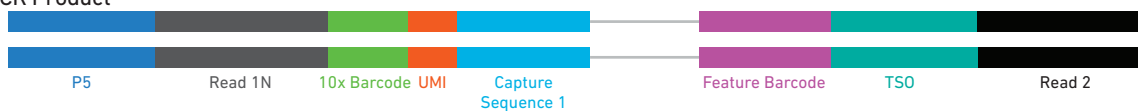

5'-AATGATACGGCGACCACCGAGATCT-ACACTCGTCGGCAGCGTCAGATGTGTATAAGAGACAG-16N-12N-TTGCTAGGACCGCCCTAAAGC-sgRNA_Feature_Barcode-CCCATGTAAGTCTGCGTTGATACCACTGCTT-AGATCGGAAGAGCACACGTCTGAACTCCAGTCAC-3'
3'-TTACTATGCCGCTGGTGGCTCTAGA-TGTGAGCAGCCGTCGCAGTCTACACATATTCTGTGTC-16N-12N-AACGATCCTGGCCGGAATTTTCG-sgRNA_Feature_Barcode-GGGTACATGAGACGCAACTATGGTACGAA-TCTAGCCTTCTCGTGTGCAGACTTGAGGTCAGTG-5'

Protocol Step 4.4 – Sample Index PCR (for CRISPR Screening Library Construction)

Feature SI
Primers 2
PN-2000099

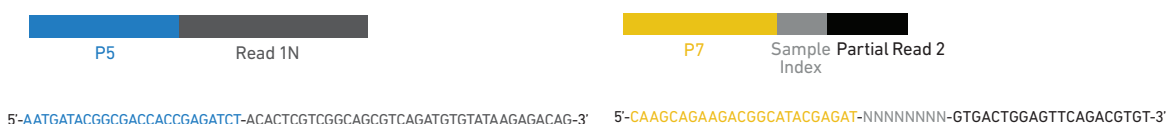

Chromium i7
Sample Index
PN-220103

Sample Index
PCR Product



5'-AATGATACGGCGACCACCGAGATCT-ACACTCGTCGGCAGCGTCAGATGTGTATAAGAGACAG-16N-12N-TTGCTAGGACCGCCCTAAAGC-sgRNA_Feature_Barcode-AGATCGGAAGAGCACACGTCTGAACTCCAGTCAC-8N-ATCTCGTATGCCGCTCTCTGCTTG-3'
3'-TTACTATGCCGCTGGTGGCTCTAGA-TGTGAGCAGCCGTCGCAGTCTACACATATTCTGTGTC-16N-12N-AACGATCCTGGCCGGAATTTTCG-sgRNA_Feature_Barcode-TCTAGCCTTCTCGTGTGCAGACTTGAGGTCAGTG-8N-TAGAGCATAACGGCAGAAGACGAAC-5'